



Peptide Test Mix

Conditions

Column: ACE 5 C18-300, 250 x 4.6mm
Part Number: ACE-221-2546
Mobile Phase: A. 0.1% TFA in H₂O
B. 0.1% TFA in MeCN
Flow Rate: 1.0ml/min
Gradient:

| T(mins) | %A | %B |
|---------|----|----|
| 0 | 90 | 10 |
| 25 | 60 | 40 |

Temperature: Ambient
Detection: UV, 220nm

Compounds

1. Gly-Tyr
2. Oxytocin
3. Angiotensin II
4. Neurotensin

