



Conditions

Columns: ACE Excel 5 C18-AR
 Dimensions: 250 x 4.6 mm
 Part Numbers: EXL-129-2546U
 Mobile Phase: MeOH/20 mM Na₂HPO₄ solution (pH 6) (15:85 v/v)
 Flow Rate: 1.0 mL/min
 Injection: 10 µL
 Temperature: Ambient
 Detection: UV, 262 nm
 Method: ChP.2015, Vol. 2


Application #AN6600

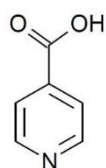


1. Impurity A

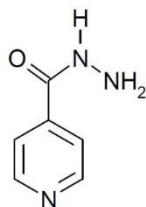


2. Isoniazid

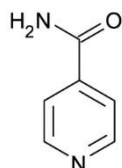


3. Impurity B

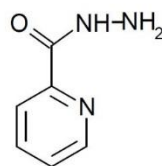

4. Impurity C

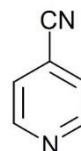

5. Impurity D

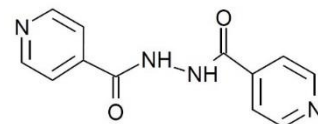

6. Impurity E

