



Angiotensin II Receptor Antagonists by LC-UV

ACE[®]
Ultra-inert
UHPLC & HPLC Columns


Application #AN3460

Conditions


Column: ACE 5 C18
Dimensions: 250 x 4.6 mm
Part Number: ACE-121-2546
Mobile Phase: 0.025 M KH_2PO_4 pH 6.0/MeCN (65:35 v/v)
Flow Rate: 1.5 mL/min
Temperature: 40 °C
Detection: UV, 220 nm

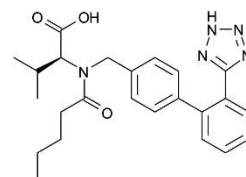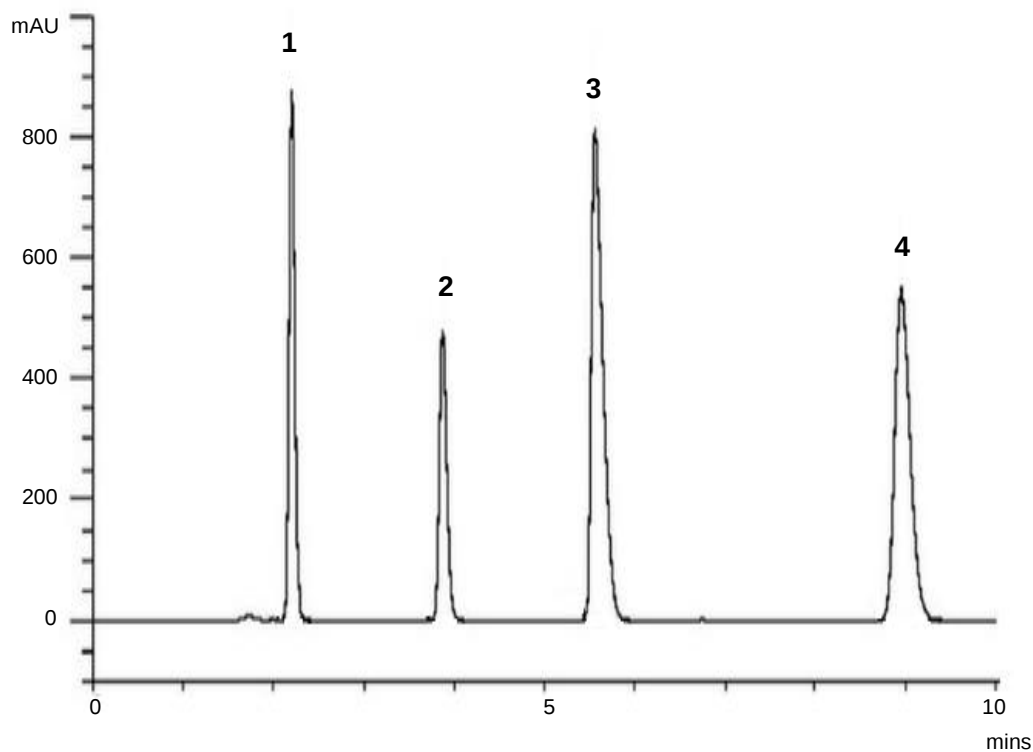


1. Valsartan

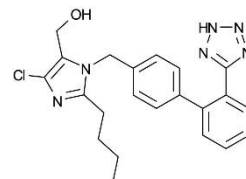

2. Losartan

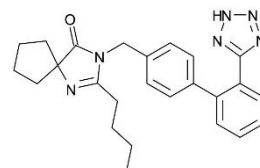

3. Irbesartan

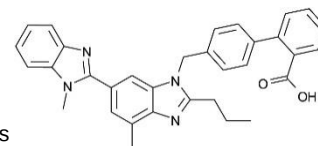

4. Telmisartan

Elshanawane AA, Abdelaziz LM, Hafez HM (2012) Stability Indicating HPLC Method for Simultaneous Determination of Several Angiotensin-II Receptor Antagonists in Their Dosage Forms. *Pharmaceut Anal Acta* 3:175. doi:10.4172/2153-2435.1000175

© 2015 ACT Ltd. All rights reserved. All products are available worldwide.

For more information contact your local ACE distributor or visit
www.ace-hplc.com or email: info@ace-hplc.com


ACE[®]
UHPLC & HPLC
Columns