


Conditions

Column: ACE UltraCore 2.5 SuperC18
 Dimensions: 100 x 3.0 mm
 Part Number: CORE-25A-1003U
 Mobile Phase: A: 20 mM ammonium formate pH 3.0 in H₂O
 B: 20 mM ammonium formate pH 3.0 in MeOH/H₂O (9:1 v/v)

Time (mins)	%B
0.0	0
7.5	100
8.5	100
9.0	0

Flow Rate: 1.2 mL/min
 Injection: 2 µL
 Temperature: 25 °C
 Detection: UV, 254 nm

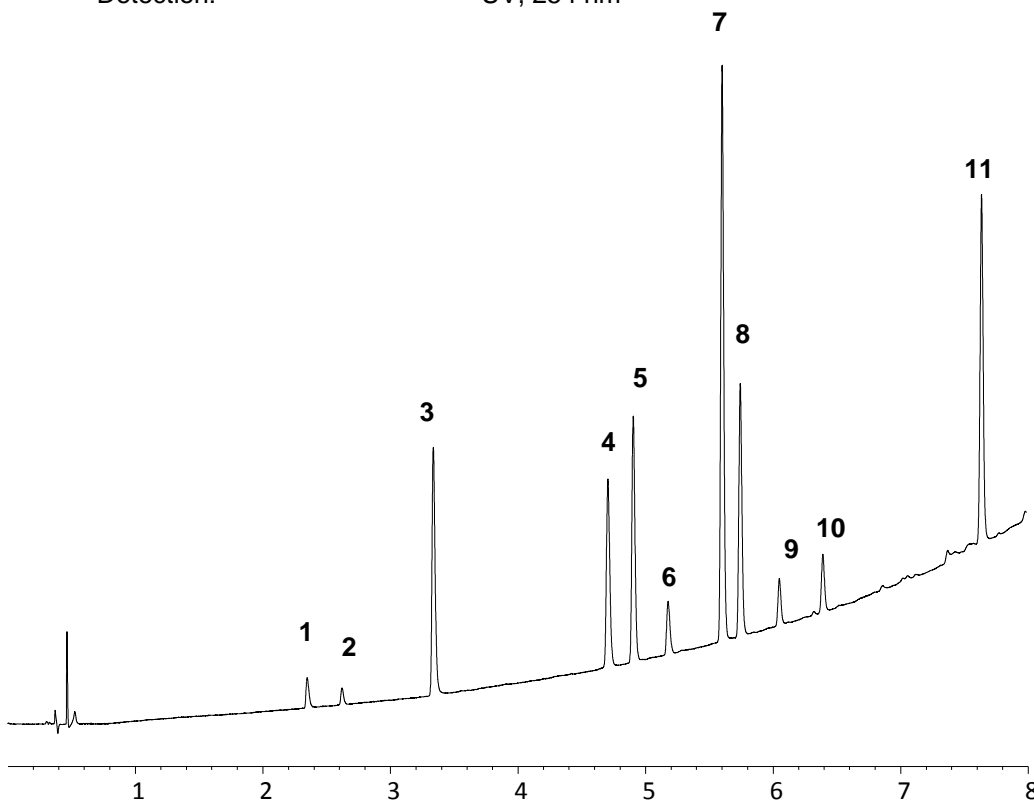

1. Pseudoephedrine
2. Scopolamine
3. Doxylamine
4. Chlorpheniramine
5. Triprolidine
6. Diphenhydramine
7. Acrivastine
8. Promethazine
9. Fexofenadine
10. Cetirizine
11. Loratadine

