


Conditions

Column: ACE Excel 1.7 CN-ES
Dimensions: 50 x 3.0 mm
Part Number: EXL-1713-0503U
Mobile Phase: A: 0.1% formic acid in H₂O
B: 0.1% formic acid in MeCN

Time (mins)	%B
0.0	5
5.0	90
6.0	90
6.5	5

Flow Rate: 0.7 mL/min
Injection: 0.5 µL
Temperature: 30 °C
Detection: UV, 240 nm

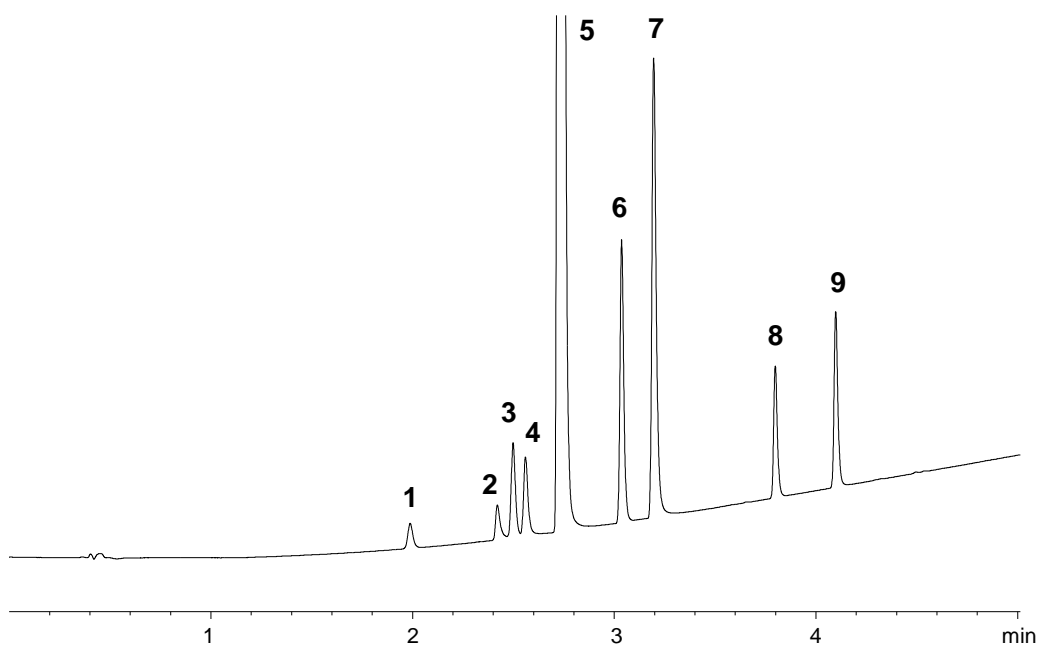

1. 2-Acetamidophenol
2. 4-Hydroxyisophthalic acid
3. Acetanilide
4. Salicylamide
5. Aspirin
6. Phenacetin
7. Salicylic acid
8. Acetylsalicylsalicylic acid
9. Salsalate

