


Corticosteroids by LC-MS/MS

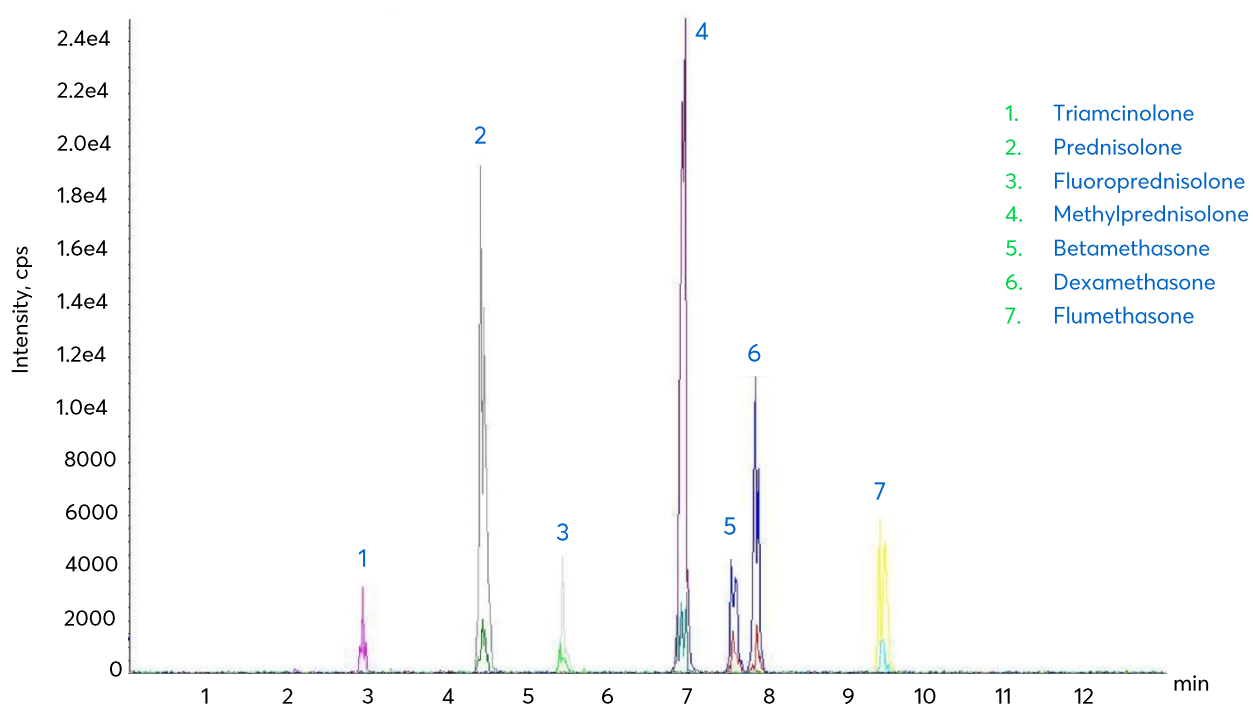

Method Details

CONDITIONS

Column:	Avantor® ACE® 3 C18-PFP										
Particle Size:	3 µm										
Dimensions:	150 x 2.1 mm										
Mobile Phase:	A: 0.1% formic acid in H ₂ O B: MeCN/0.1% formic acid in H ₂ O (90:10 v/v)										
Gradient:	<table border="1"> <thead> <tr> <th>Time (mins)</th><th>%B</th></tr> </thead> <tbody> <tr> <td>0.0</td><td>30</td></tr> <tr> <td>14.0</td><td>50</td></tr> <tr> <td>17.0</td><td>95</td></tr> <tr> <td>20.0</td><td>30</td></tr> </tbody> </table>	Time (mins)	%B	0.0	30	14.0	50	17.0	95	20.0	30
Time (mins)	%B										
0.0	30										
14.0	50										
17.0	95										
20.0	30										
Flow Rate:	0.3 mL/min										
Injection:	25 µL										
Temperature:	15 °C										
Detection:	Turbospray, MRM										

ORDERING TABLE

Product	Details	Size	Part number
Avantor® ACE® 3 C18-PFP	HPLC Column	150 x 2.1 mm	ACE-1110-1502

Reproduced with permission of Laboratorio Salud Publica de Burgos, Spain