



Cyclodextrin-Encapsulated Flavour Compounds in Beer

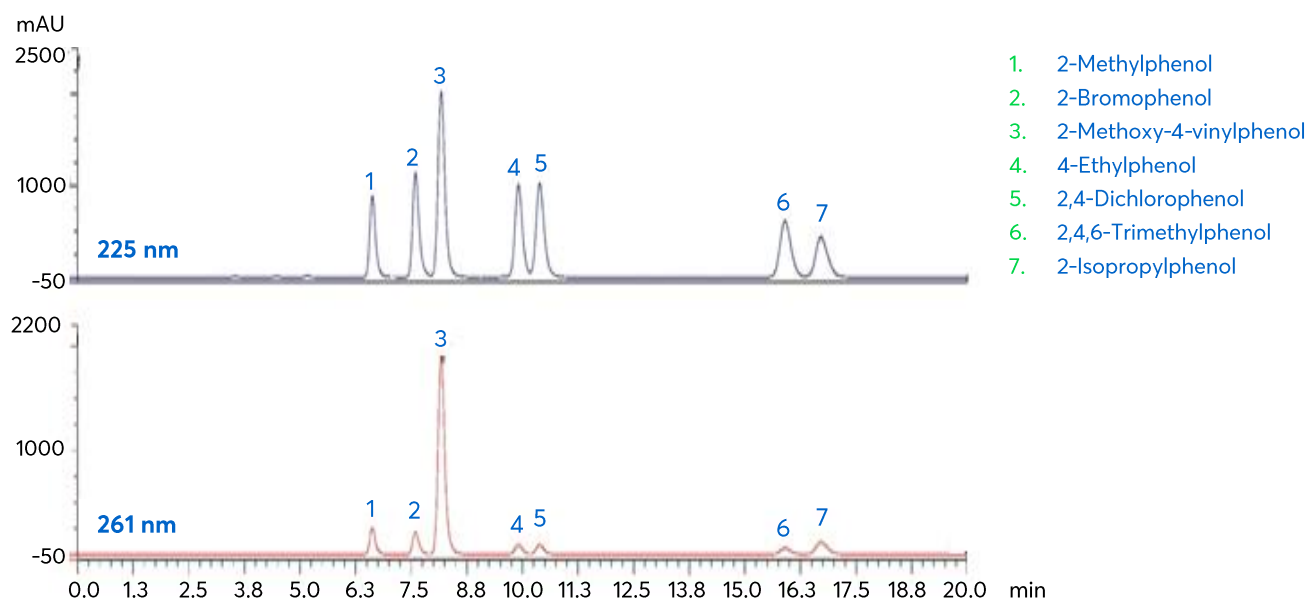

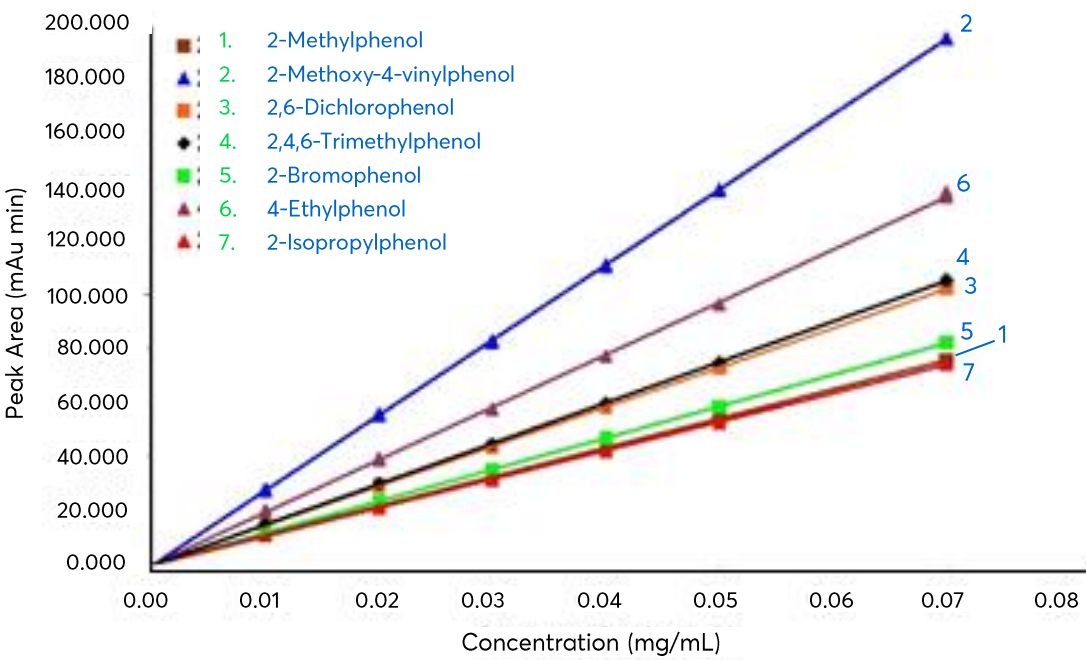

Method Details

CONDITIONS

Column:	Avantor® ACE® 3 C18
Particle Size:	3 µm
Dimensions:	150 x 4.0 mm
Mobile Phase:	0.1% phosphoric acid in MeOH/H ₂ O (53:47 v/v)
Flow Rate:	0.5 mL/min
Injection:	20 µL
Temperature:	35 °C
Detection:	UV, 225 nm (261 nm for 2-Methoxy-4-vinylphenol)

ORDERING TABLE

Product	Details	Size	Part Number
Avantor® ACE® 3 C18	HPLC Column	150 x 4.0 mm	ACE-111-1504

Reproduced with permission of FlavorActiV Ltd, UK