


Determination of Baicalin in Children's Liquid Oral Medicine

ACE[®]
Ultra-inert
UHPLC & HPLC Columns


Application #AN6290

Conditions

Column: ACE Excel 5 AQ
Dimensions: 150 x 4.6 mm
Part Number: EXL-126-1546U
Mobile Phase: 0.1% H₃PO₄ in H₂O/ MeOH (50:50 v/v)
Flow Rate: 1.0 mL/min
Injection: 5 µL
Temperature: 40 °C
Detection: UV, 280 nm
Monograph Method: ChP.2015, Vol. 1

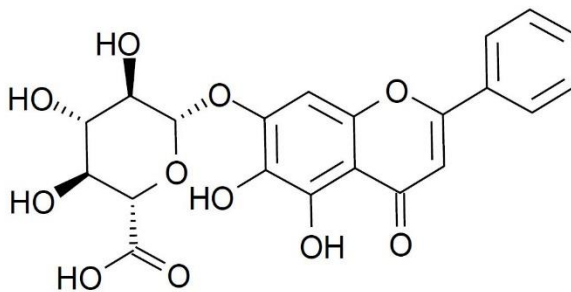

1. Baicalin

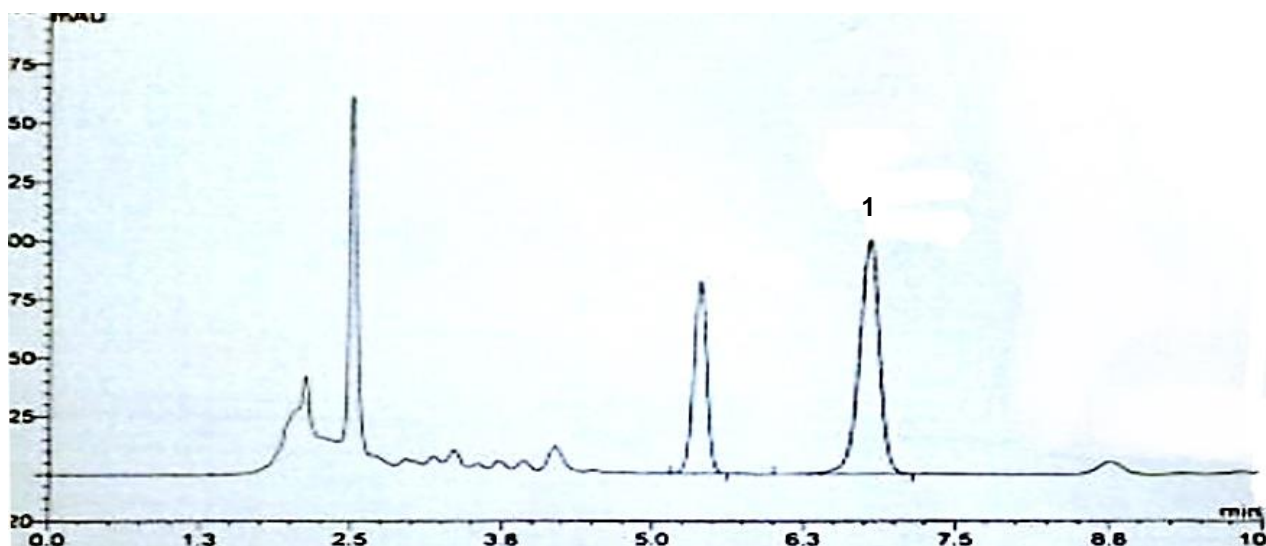

© 2018 ACT Ltd. All rights reserved. All products are available worldwide.

For more information contact your local ACE distributor or visit
www.ace-hplc.com or email: info@ace-hplc.com


ACE[®]
UHPLC & HPLC
Columns