






Digitoxin

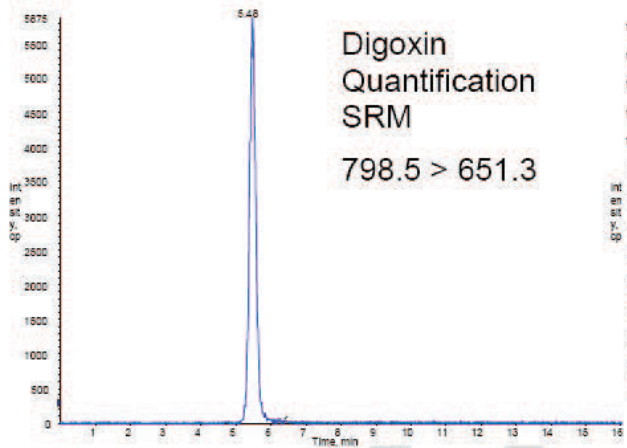


Digitoxin
Quantification
SRM
798.5 > 651.3

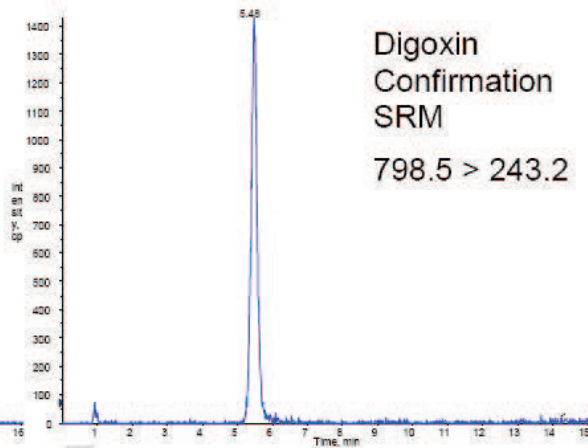

Digitoxin
Confirmation
SRM
798.5 > 243.2

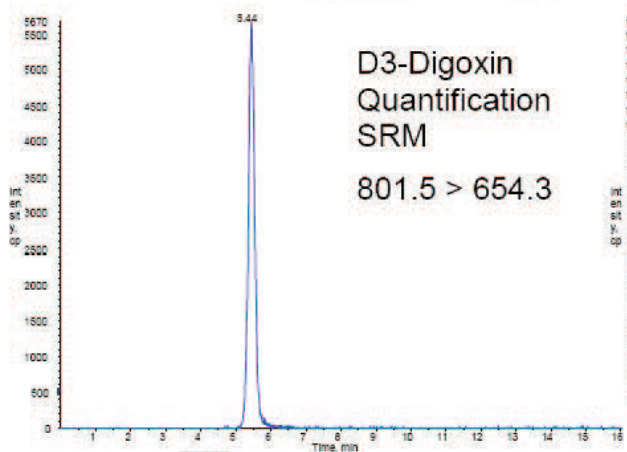

D3-Digoxin
Quantification
SRM
801.5 > 654.3

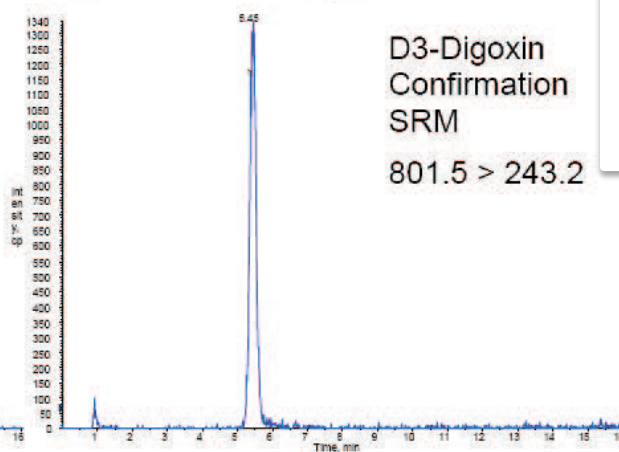

D3-Digoxin
Confirmation
SRM
801.5 > 243.2

Fortis C18 100x2.1mm 3um

0.1% formic acid in Water
0.1% formic acid in MeOH

30:70 0mins
70:30 10mins

0.2ml/min

30oC

MS

Digitoxin
D3-Digitoxin