
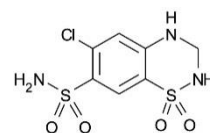


Conditions


Column: ACE Excel 2 C18-PFP
 Dimensions: 50 x 3.0 mm
 Part Number: EXL-1010-0503U
 Mobile Phase: A: 10 mM ammonium formate pH 3.0 in H₂O
 B: 10 mM ammonium formate pH 3.0 in MeOH/H₂O (9:1 v/v)

Time (mins)	%B
0.0	5
0.5	5
5.0	70
5.5	70
6.0	5


Flow Rate: 1 mL/min
 Injection: 2 µL
 Temperature: 60 °C
 Detection: UV, 254 nm

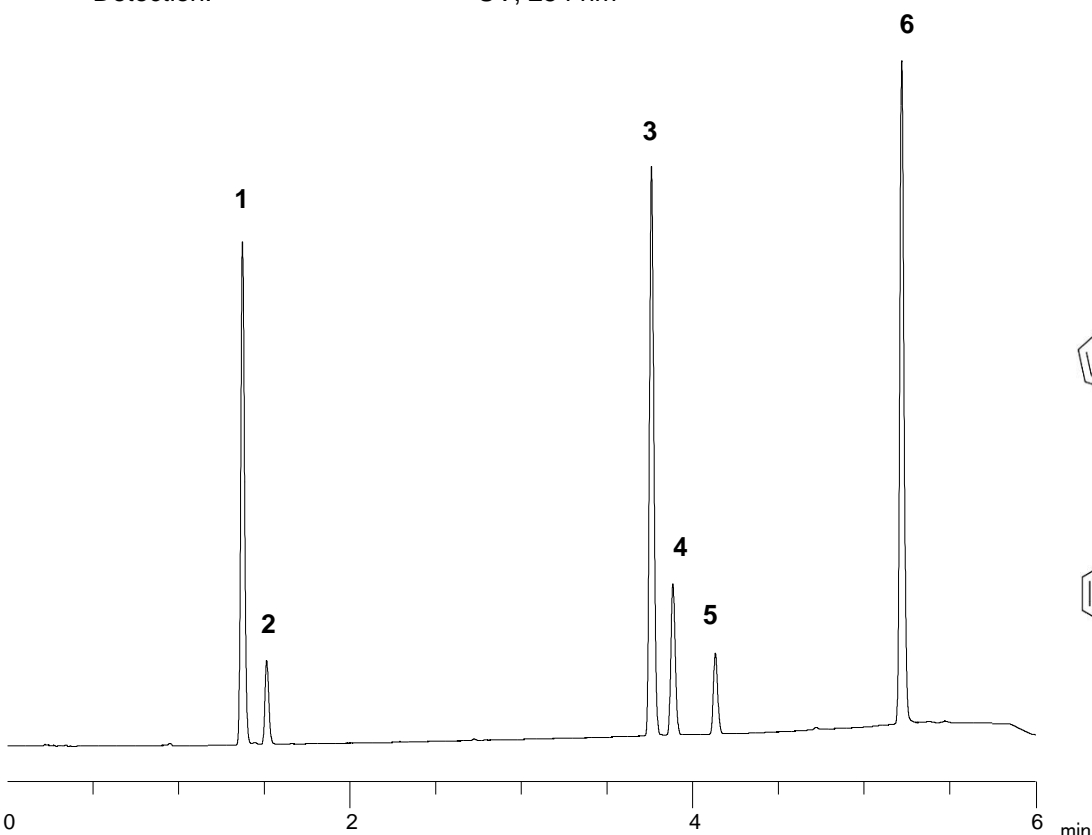


1. Hydrochlorothiazide

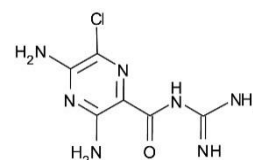


2. Amiloride

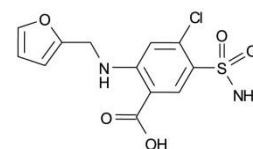


3. Furosemide

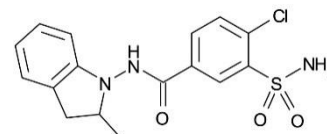

4. Indapamide

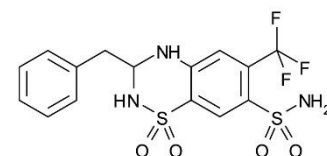

5. Bendroflumethiazide

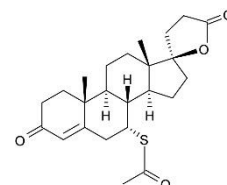

6. Spironolactone