

Exploring The Power Of Chromatographic Selectivity For Polar & Non-Polar Analytes With A Unique HPLC / UHPLC Polar Embedded Stationary Phase


Alan P McKeown¹, Geoffrey Faden²

¹Advanced Chromatography Technologies Ltd, 1 Berry Street, Aberdeen, Scotland, AB25 1HF UK ²MACMOD Analytical Inc., 103 Commons Court, PO Box 587, Chadds Ford, PA 19317 USA


THE CHALLENGE

Engineer a new, unique HPLC / UHPLC phase with polar retention / resolution and alternative selectivity that is reproducible, robust and gives efficient chromatography

1. RESOLUTION, SELECTIVITY, EFFICIENCY & RETENTION

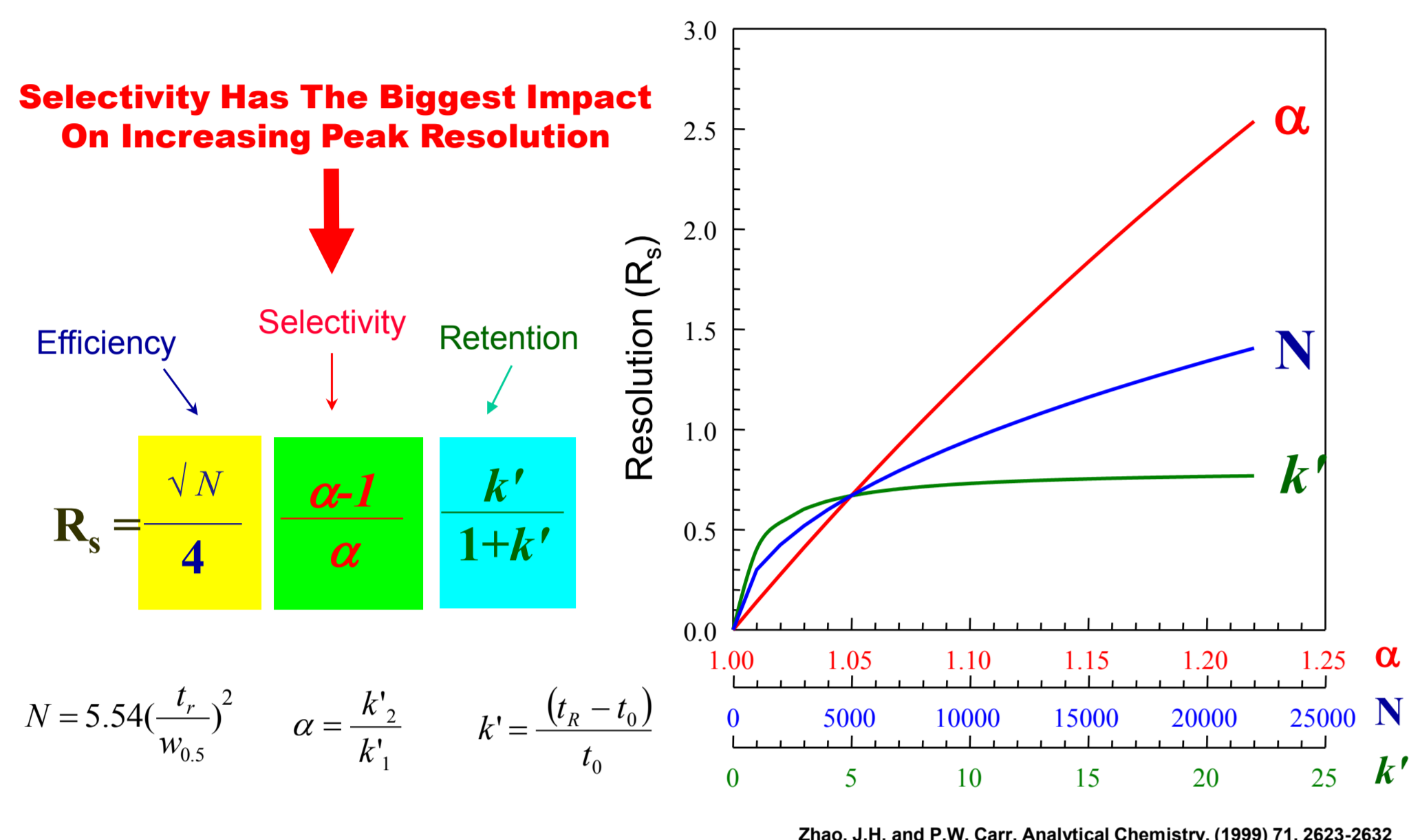


2. C18 STATIONARY PHASES SHOW SIMILAR SELECTIVITY

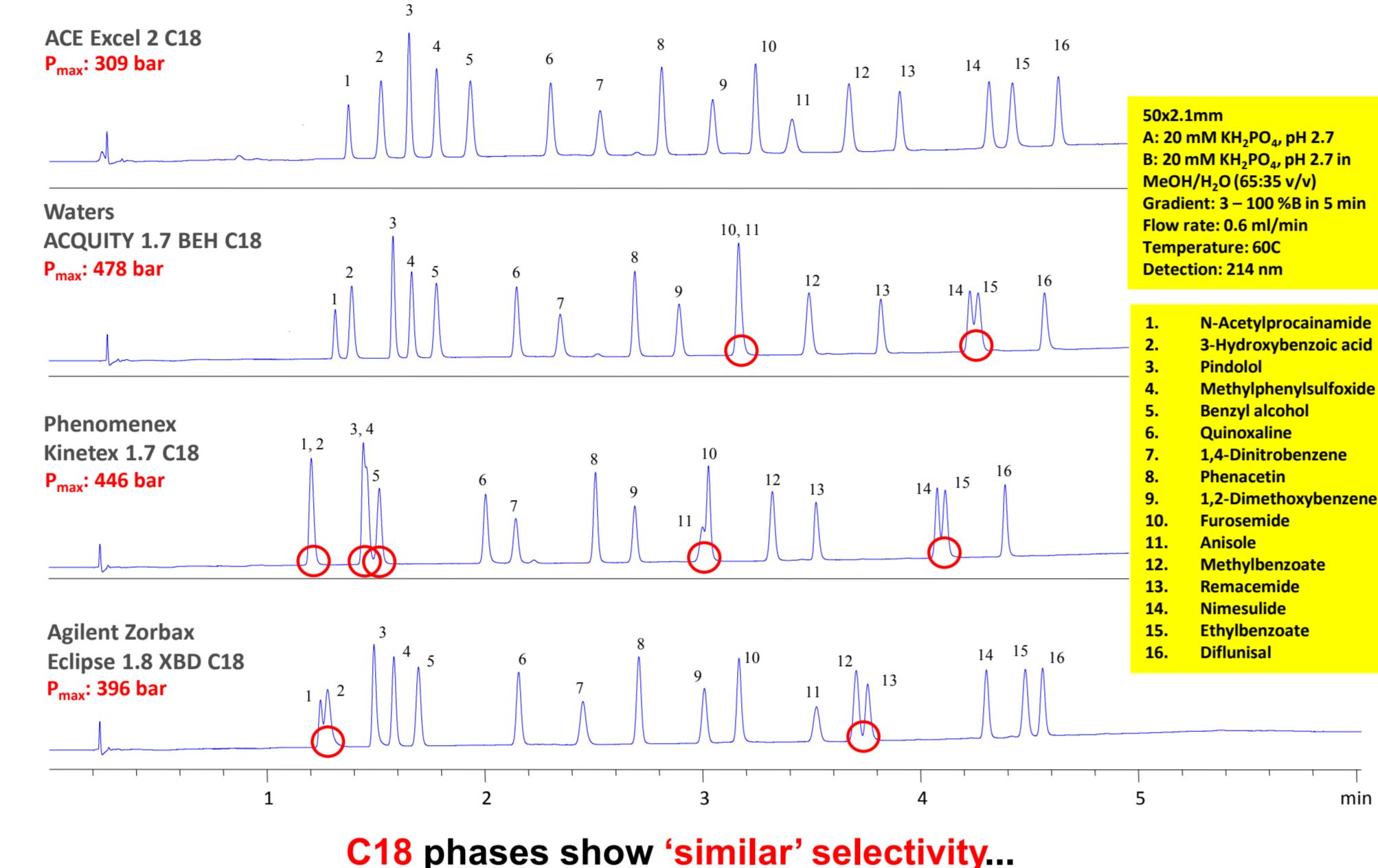

3. ACE® C18-Amide™: A NEW POLAR EMBEDDED PHASE OPTION


- Uniquely designed ligand (USP L1 / L60)
 - Maximise stability and multiple modes of interaction
- Enhanced retention / resolution of polar acidic analytes
 - Ideal for H-bond donor analytes: acids, amino, amides etc
- Enhanced retention and resolution for phenolics
 - Wine acids, green teas, hydroxylated / polar analytes etc
- Usable in 100% aqueous eluents: no dewetting

4. DETERMINING SELECTIVITY VALUES* FOR PHASES

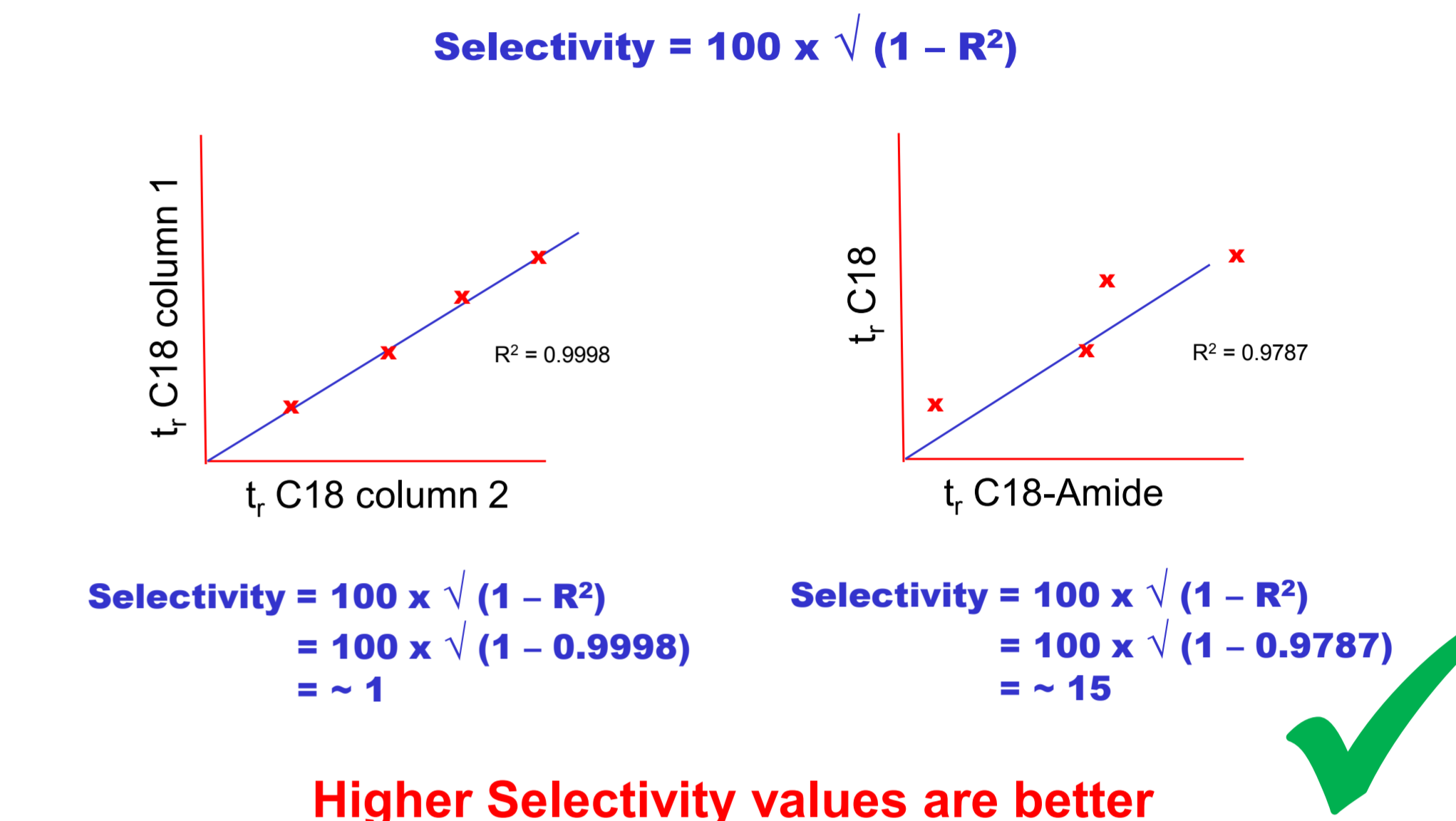


* Neue, O'Gara, Méndez "Selectivity in Reversed-Phase Separations: Influence of the Stationary Phase", J. Chromatogr. A 1127 (2006), 161-174

5. ACE® C18-Amide™ EXCELLENT POLAR SELECTIVITY

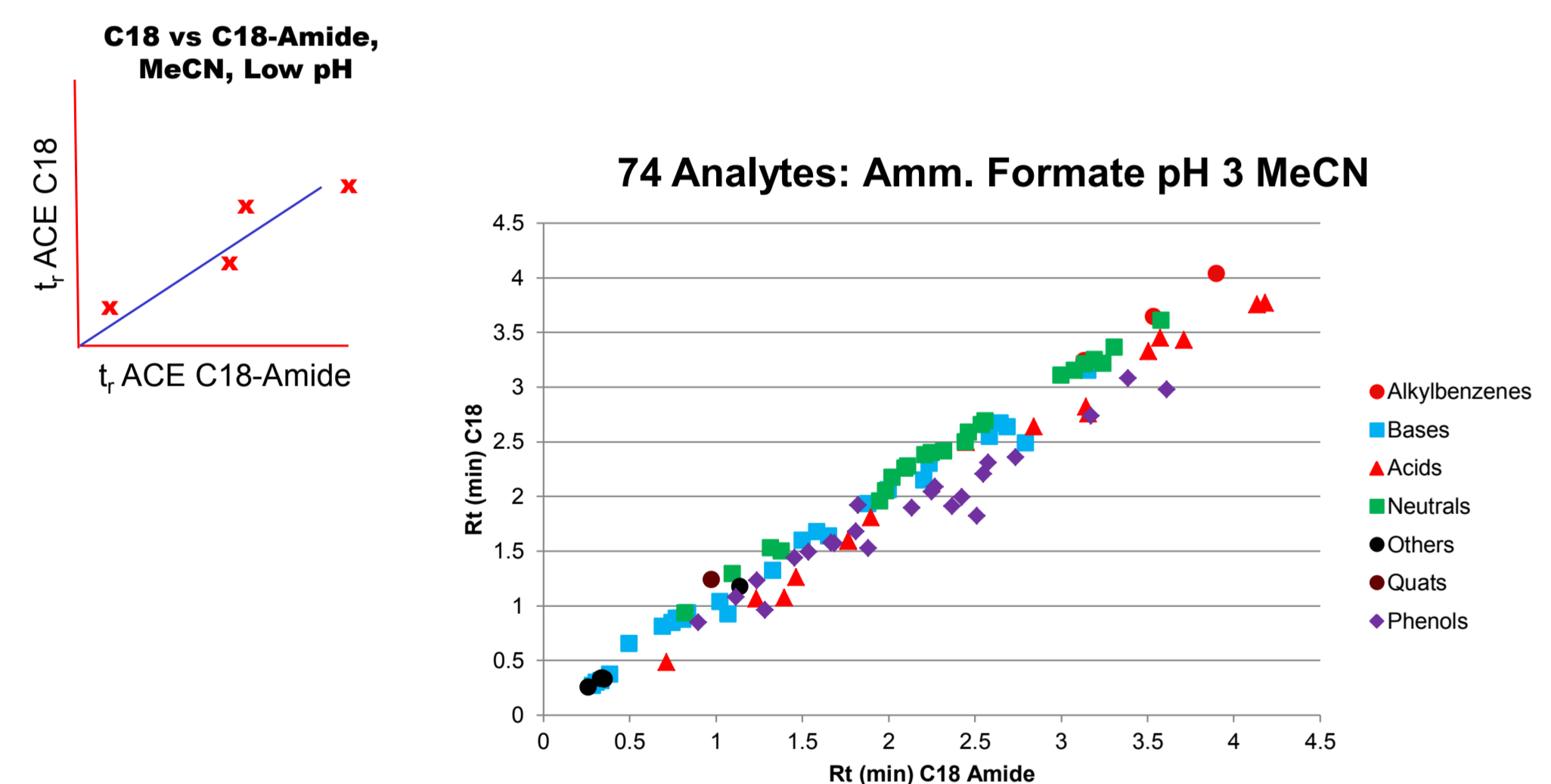

Selectivity = 20 → Powerful For Method Development


6. POLAR SELECTIVITY INCLUDING LOW & MID pH EFFECTS

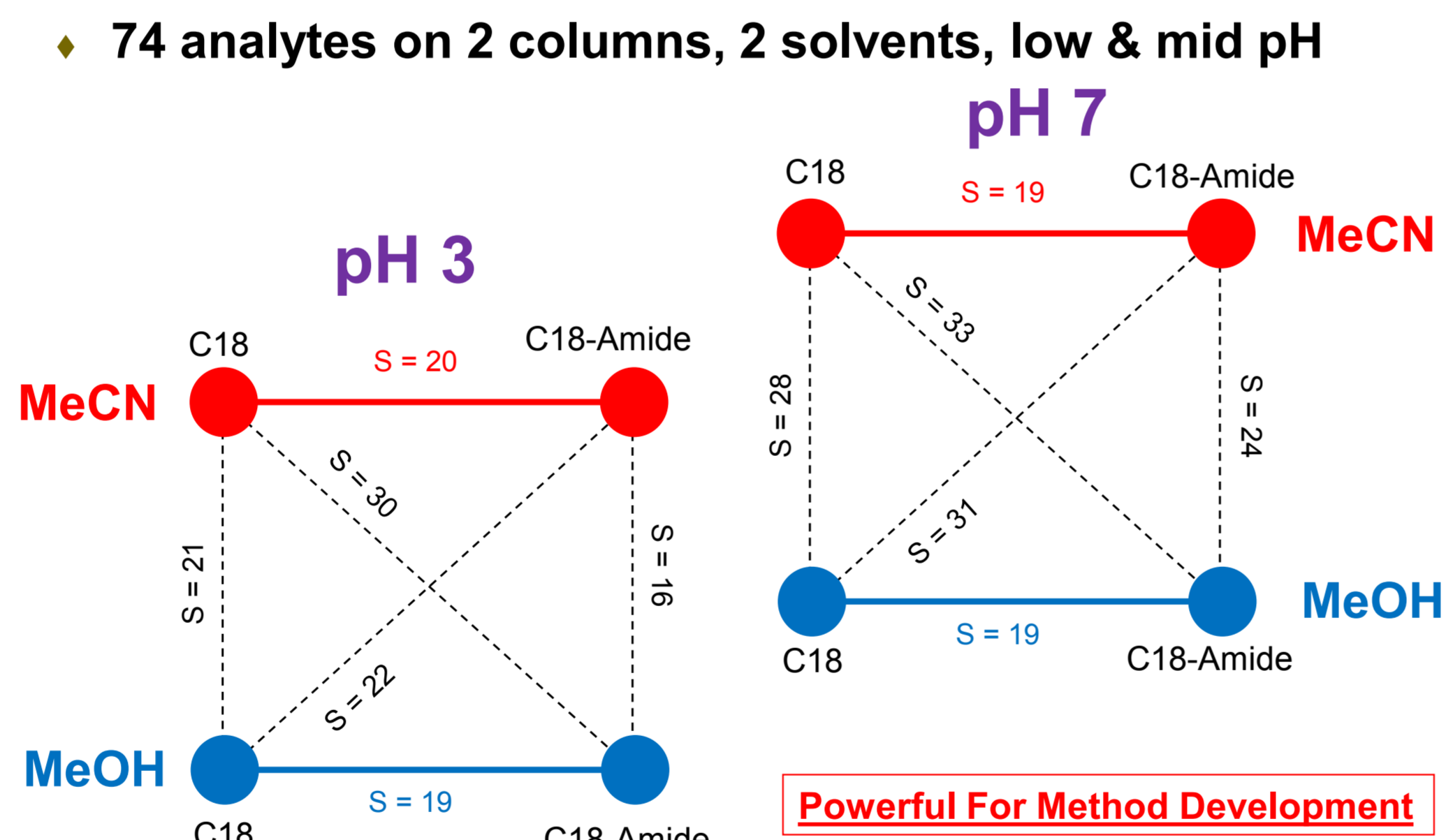


7. EMBEDDED PHASE STABILITY AT LOW & MID pH

- Real time stability data collected on retention and efficiency (acidic, neutral, basic analytes):
 - 20,000 column volumes @ 60C / pH 2.5 (phosphate)
 - 20,000 column volumes @ 60C / pH 7.0 (phosphate)


Equates to > 60 days of use (8hrs use per day)

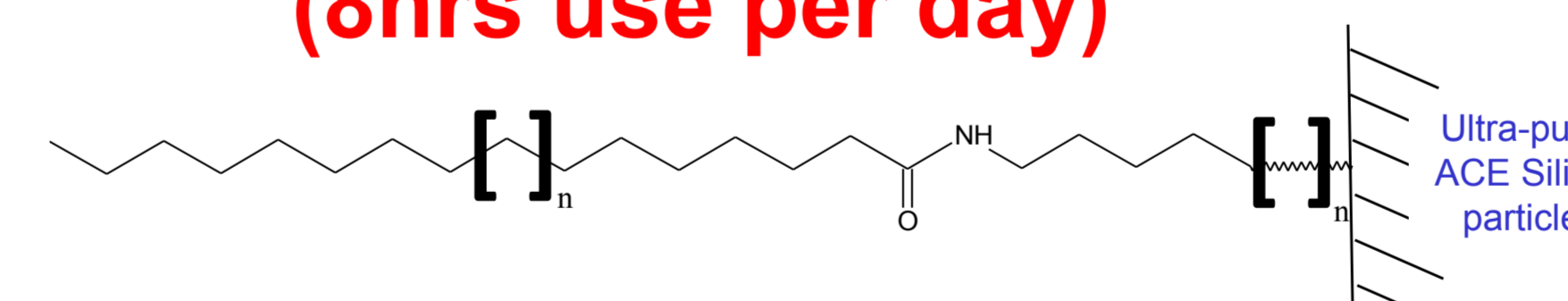


8. METHOD DEVELOPMENT: ALTERNATE POLAR SELECTIVITY

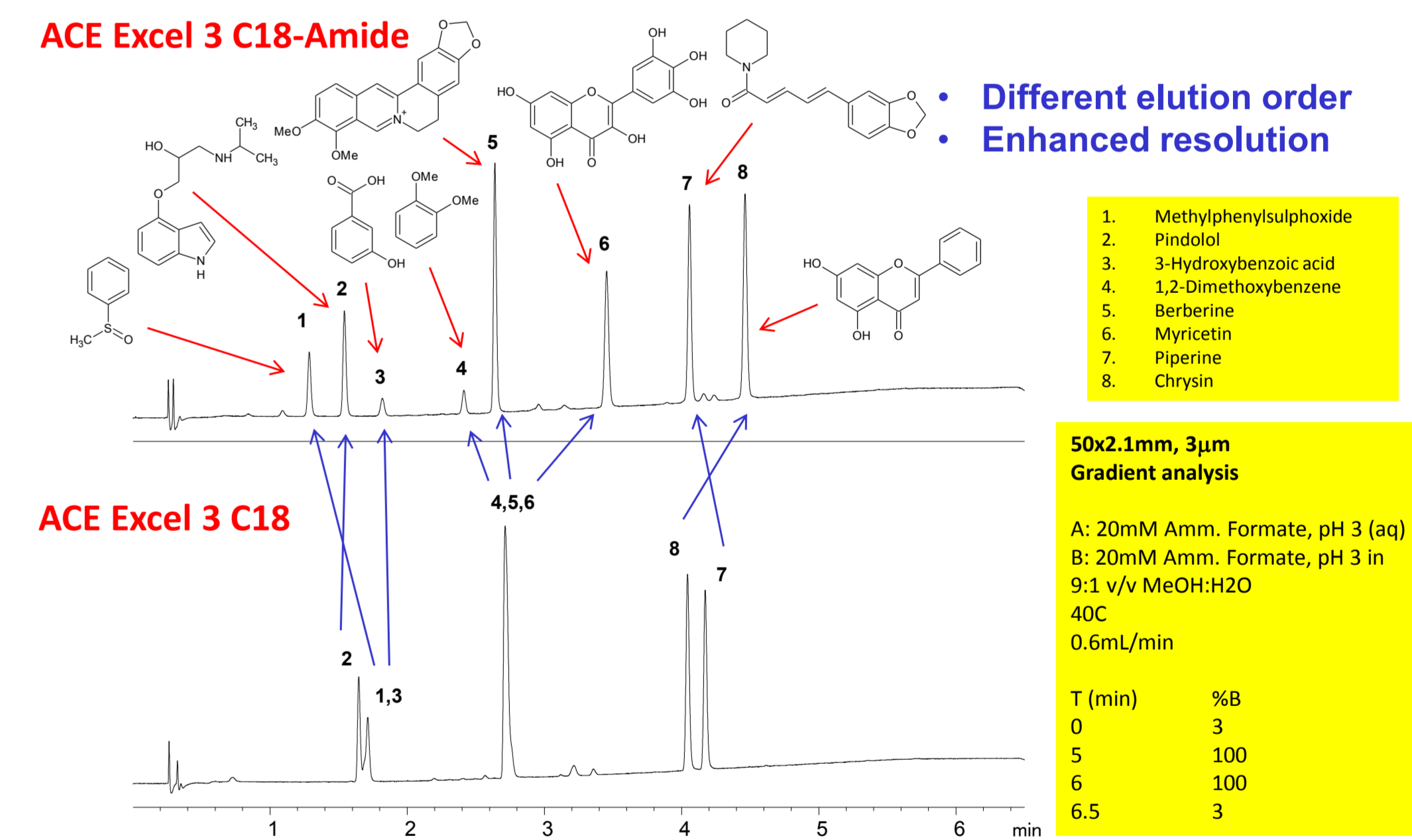


9. SEPARATION OF CATECHOLS AND RESORCINOLS

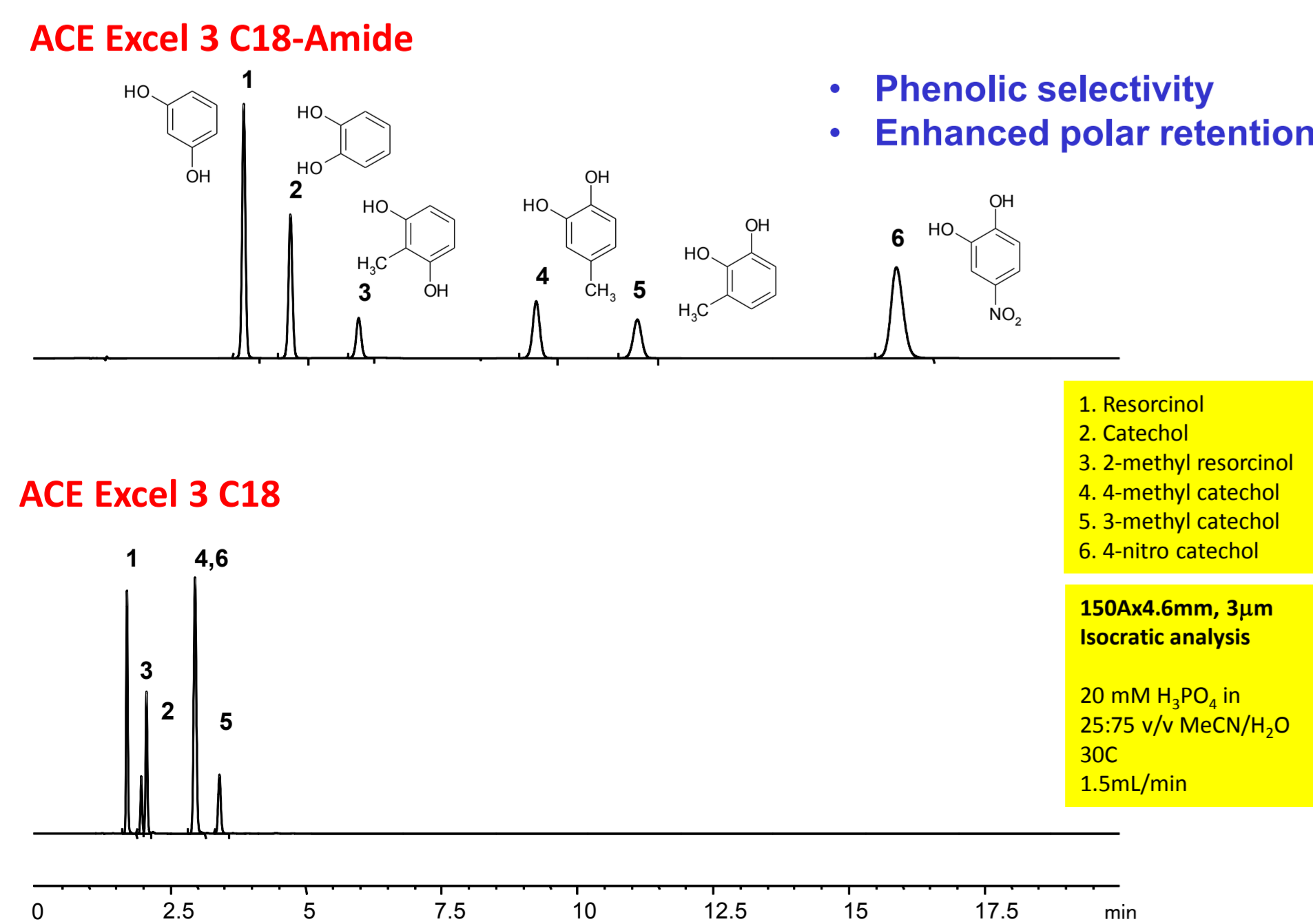

10. RAPID BEVERAGE ANALYSIS: WINE ACIDS – 100% AQUEOUS

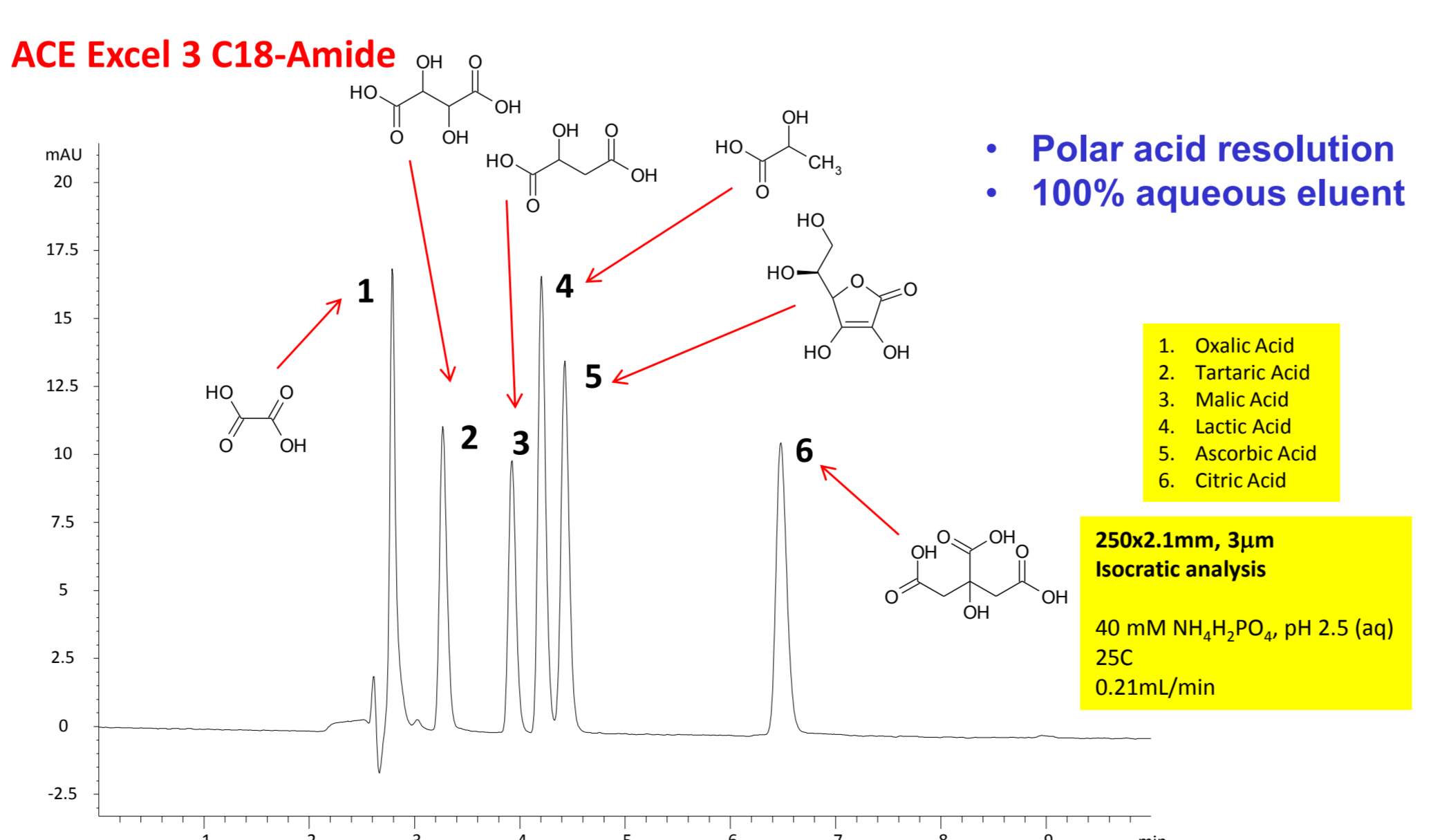

11. PHARMA COPEIPEL RELATED ANALYSES: STEROIDS

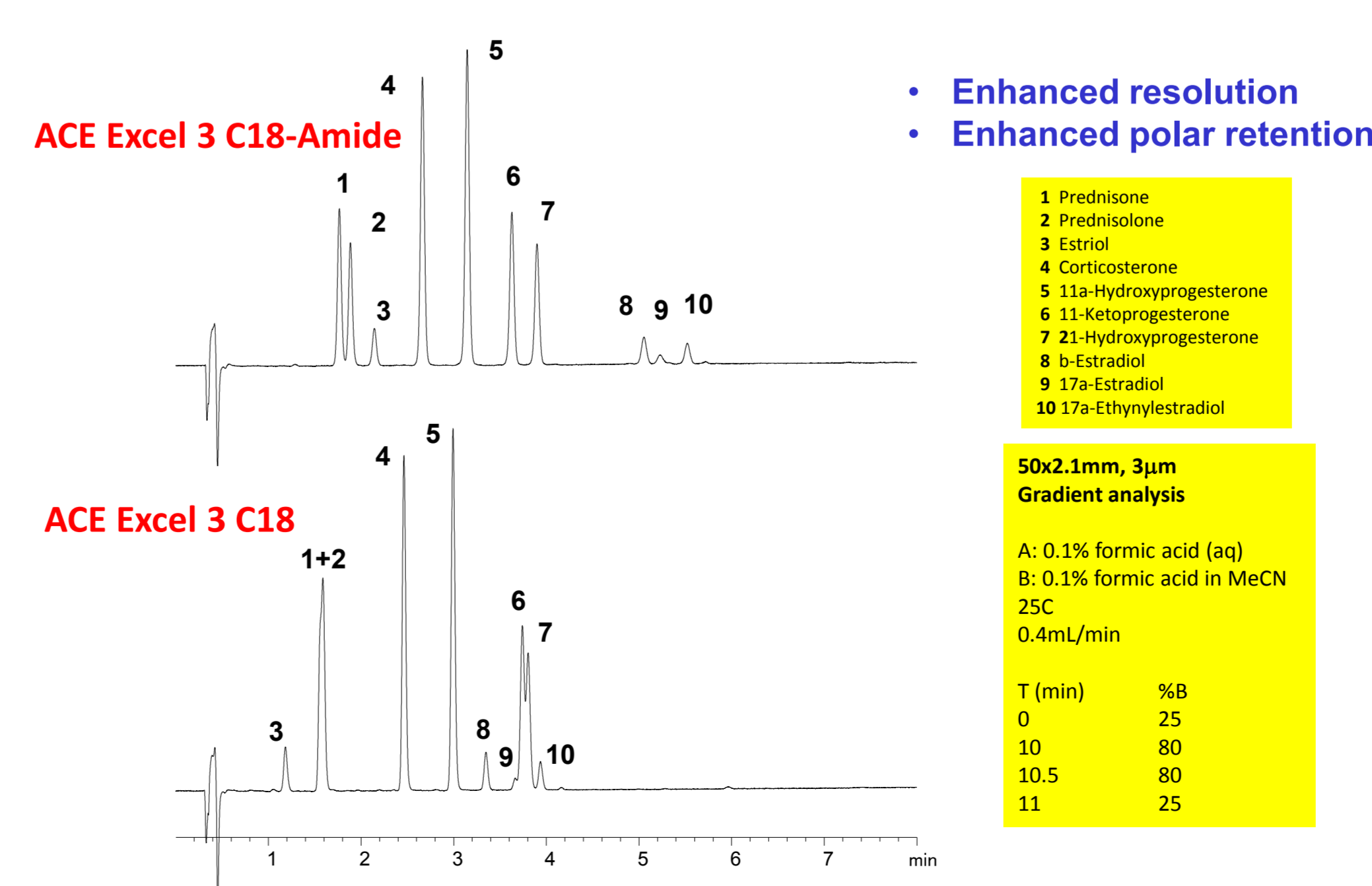

12. SUMMARY AND CONCLUSIONS

- Separations of very polar to non polar species are achievable using the ACE® C18-Amide™.
- The ACE® C18-Amide™ provides alternative selectivity to C18 based phases which is ideal for method development or sample screening purposes. Further applications are available.
- The unique ligand design of the ACE® C18-Amide™ improves the hydrophobic retention mechanism contribution to separations whilst providing enhanced stationary phase stability.
- The ACE® C18-Amide™ provides chromatographers and method developers with a NEW selectivity option for mixtures containing very polar and / or non-polar analytes.

ACEB is a registered trademark of Advanced Chromatography Technologies Limited. ACE Excel™ is a trademark of Advanced Chromatography Technologies Limited. Advanced Chromatography Technologies Limited acknowledges the registered and unregistered trademarks of Agilent Technologies Inc., Phenomenex Inc. and Waters Corporation and has no affiliation with any of these companies.