


Garlic Analysis (II)

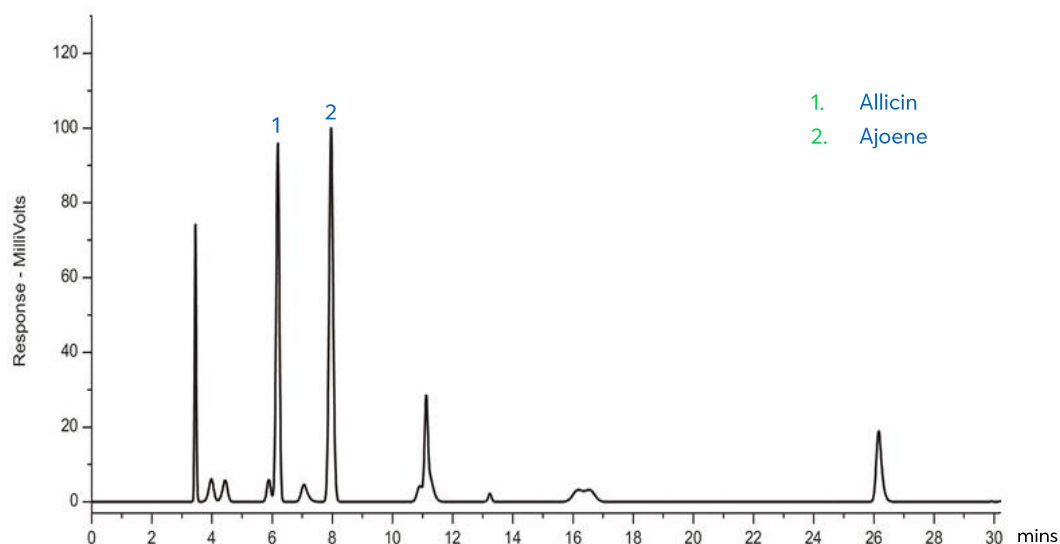

Method Details

CONDITIONS

Column: Avantor® ACE® 5 C18

Particle Size: 5 µm

Dimensions: 250 x 4.6 mm

Mobile Phase: A: H₂O
B: MeCN

Gradient:


Time (mins)	% B
0	40
20	100
25	100

Flow Rate: 1 mL/min


Injection: 20 µL

Temperature: 30 °C

Detection: UV, 254 nm

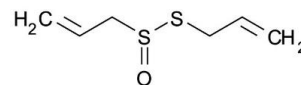

Allicin

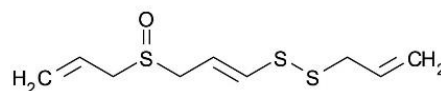

Ajoene

ORDERING TABLE

Product	Details	Size	Part Number
Avantor® ACE® 5 C18	HPLC Column	250 x 4.6 mm	ACE-121-2546

Reproduced with permission of Neem Biotech, Cardiff, UK