

Glitoxin from *Aspergillus Fumigatus* Liquid Culture

ACE[®]
Ultra-inert
UHPLC & HPLC Columns



Application #AN3780

Conditions

Column: ACE 5 C18
Dimensions: 250 x 4.6 mm
Part Number: ACE-121-2546
Mobile Phase: A: 0.05% TFA in H₂O/MeCN (90:10 v/v)
B: 0.05% TFA in H₂O/MeCN (40:60 v/v)

Time (mins)	%B
0	10
21	100

Flow Rate: 1 mL/min
Detection: UV, 254 nm

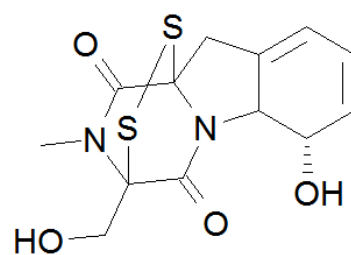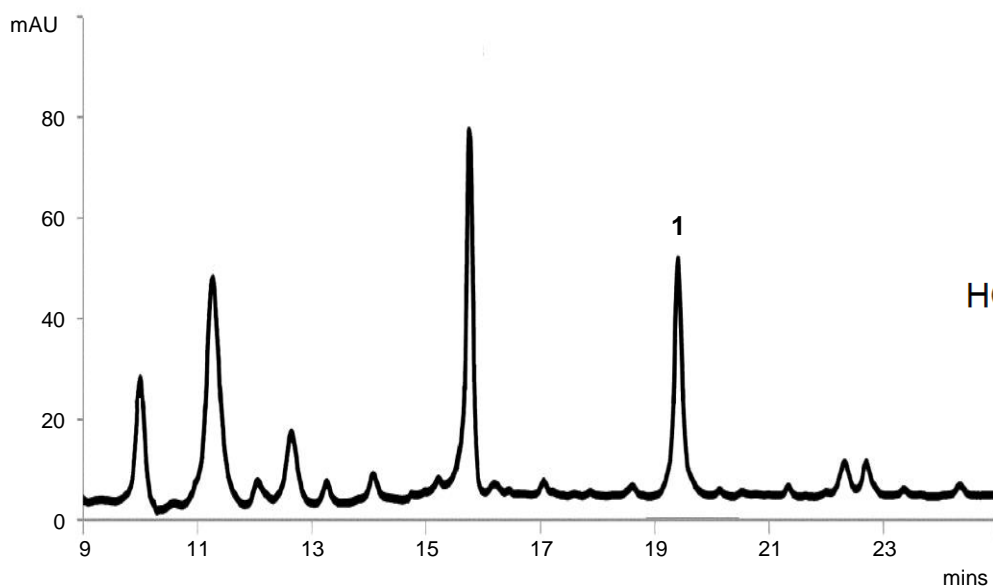

1. Glitoxin

Svahn KS, Goransson U, Chryssanthou E, Olsen B, Sjolín J, Stromstedt A. Induction of Glitoxin Secretion in *Aspergillus fumigatus* by Bacteria-Associated Molecules. PloS ONE 9(4): e93685.
doi:10.1371/journal.pone.0093685

© 2015 ACT Ltd. All rights reserved. All products are available worldwide.

For more information contact your local ACE distributor or visit
www.ace-hplc.com or email: info@ace-hplc.com

ACE[®]
UHPLC & HPLC
Columns