
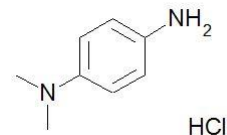



Application #AN2100

Conditions


Column: ACE Excel 5 C18-Amide
Dimensions: 250 x 4.6 mm
Part Number: EXL-1212-2546U
Mobile Phase: 1.8 g disodium phosphate dodecahydrate + 2.8 g potassium dihydrogen phosphate + 1.0 g sodium 1-heptanesulfonate (all diluted to 1.0 L with water): MeCN (60:40 v/v)
Flow Rate: 1 mL/min
Temperature: 60 °C
Detection: UV, 280 nm

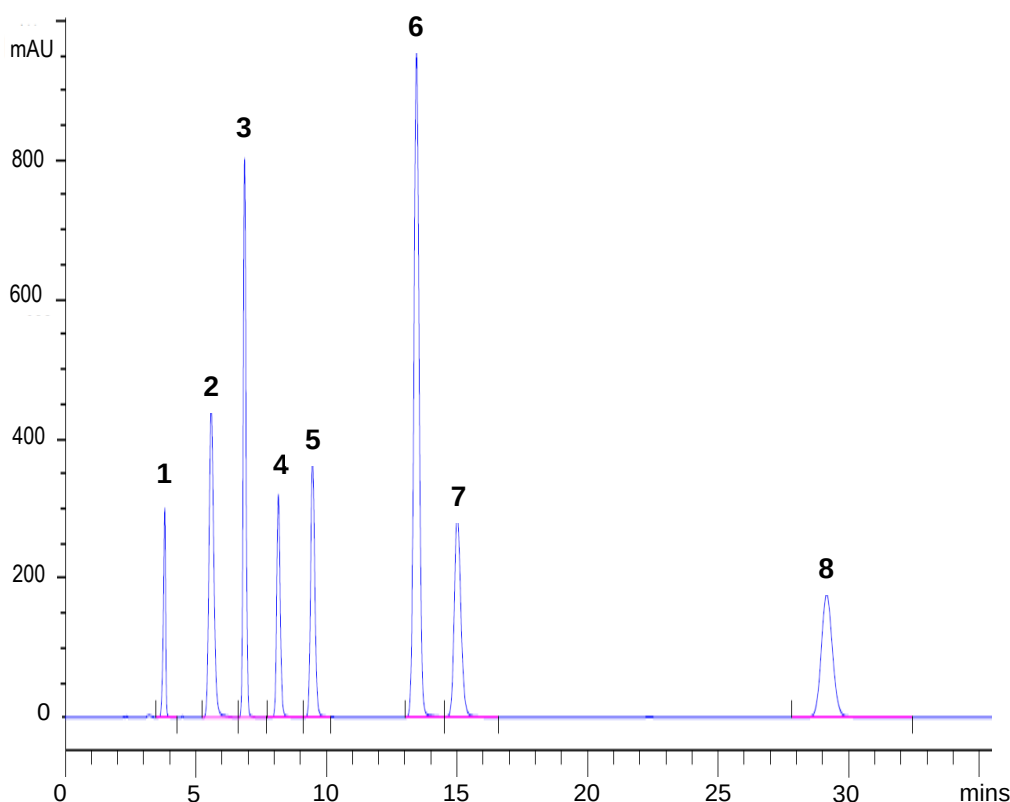


1. N,N-Dimethyl-1,4-phenylenediamine HCl

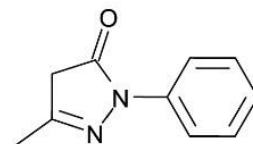


2. 5-Methyl-2-phenyl-1,2-dihydropyrazol-3-one

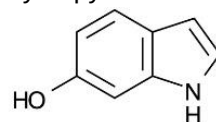


3. 6-Hydroxyindole

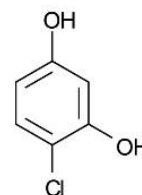


4. 4-Chlororesorcinol

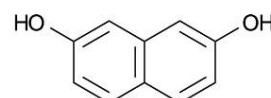


5. 2,7-Dihydroxynaphthalene

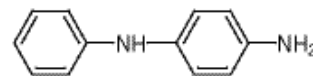

6. 4-Aminodiphenylamine

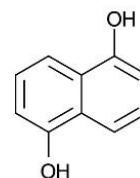

7. 1,5-Dihydroxynaphthalene

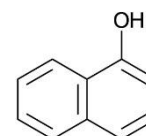

8. 1-Naphthol