

# Highly Sensitive and Selective Quantification of Microcystin Toxins in Drinking Water by UHPLC-MS/MS

Alan P McKeown<sup>1</sup>, Geoffrey Faden<sup>2</sup>,

<sup>1</sup>Advanced Chromatography Technologies Ltd, 1 Berry Street, Aberdeen, Scotland, AB25 1HF UK <sup>2</sup>MACMOD Analytical Inc., 103 Commons Court, PO Box 587, Chadds Ford, PA 19317 USA


## 1. BACKGROUND

- Cyanobacteria / blue-green algae are common. Microcystins are cyclic heptapeptides produced by cyanobacteria and associated with human hepatotoxicity.
- Under nutrient-rich conditions, cyanobacteria can rapidly accumulate and form blooms.
- Due to their hepatotoxic nature, the WHO has established a provisional limit of 1µg/L in water supplies for total Microcystin-LR as a marker for cyanobacteria toxin levels.
- This poster describes a method for low ppb quantification of Microcystin-LR, Microcystin-RR and Microcystin-YR by UHPLC-MS/MS using an ACE Excel 2µm C18 column.


## 2. ALGAE BLOOMS AND MICROCYSTINS



[http://www.aquarius-systems.com/Entries/View/349/bluegreen\\_algae.aspx](http://www.aquarius-systems.com/Entries/View/349/bluegreen_algae.aspx)

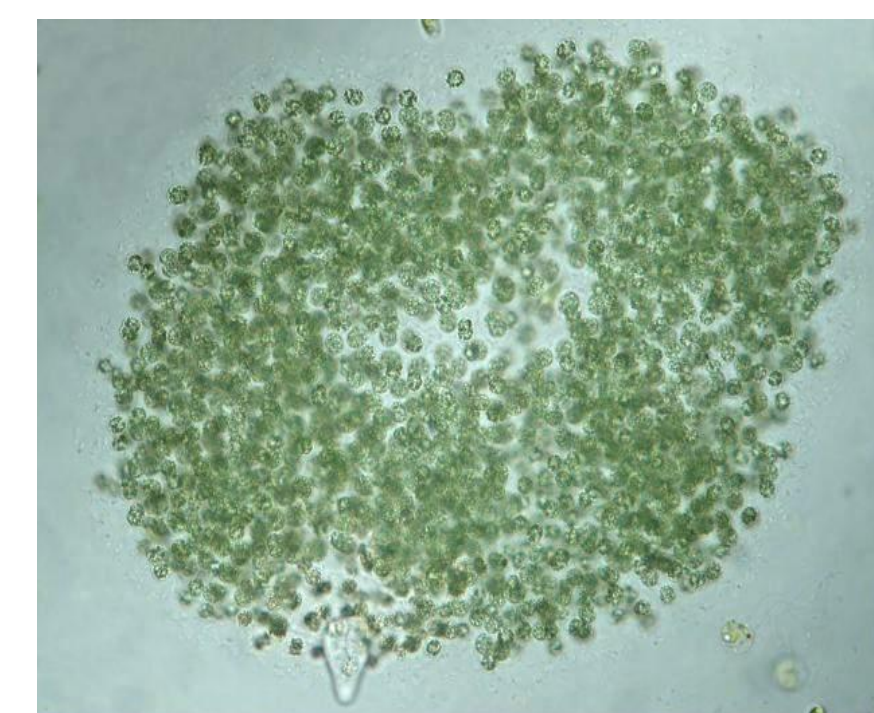


<http://www.pimgfactory.de/index.html>

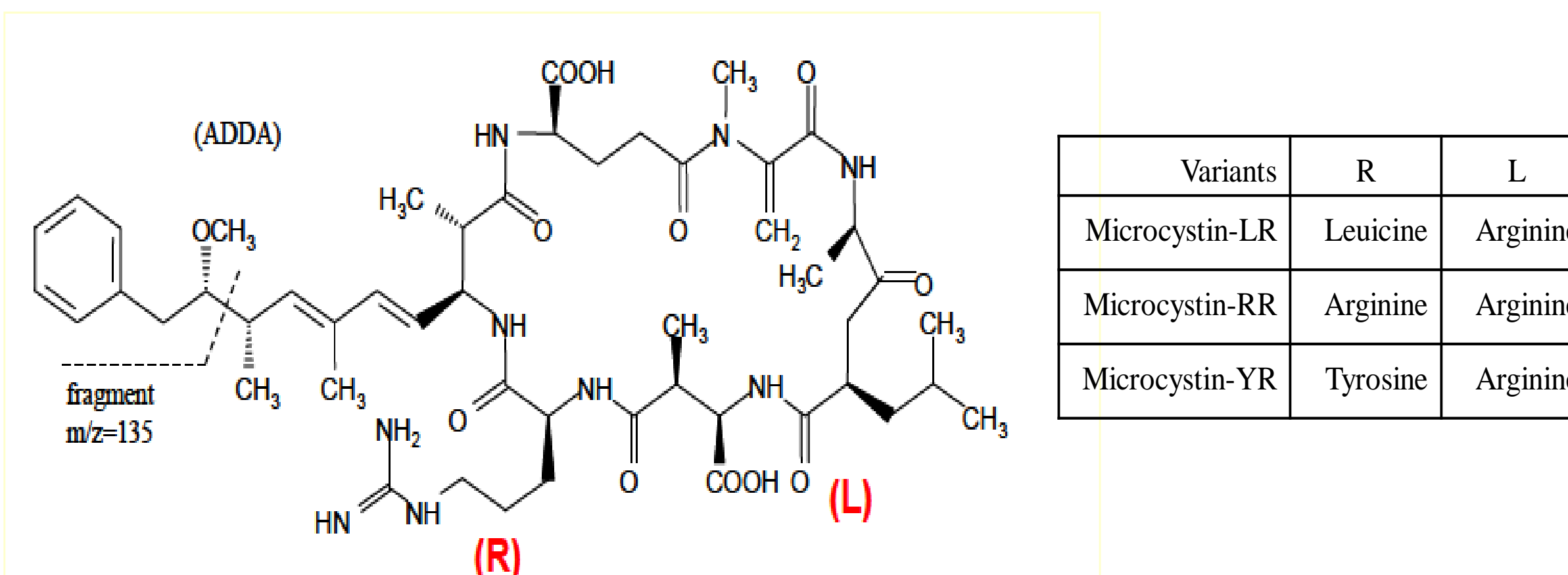


Over 80 known microcystin variants have been reported

Int. J. Environ. Res. Public Health 2012, 9

## 4. MS SCANS AND SENSITIVITY INVESTIGATIONS AT 100 PPB

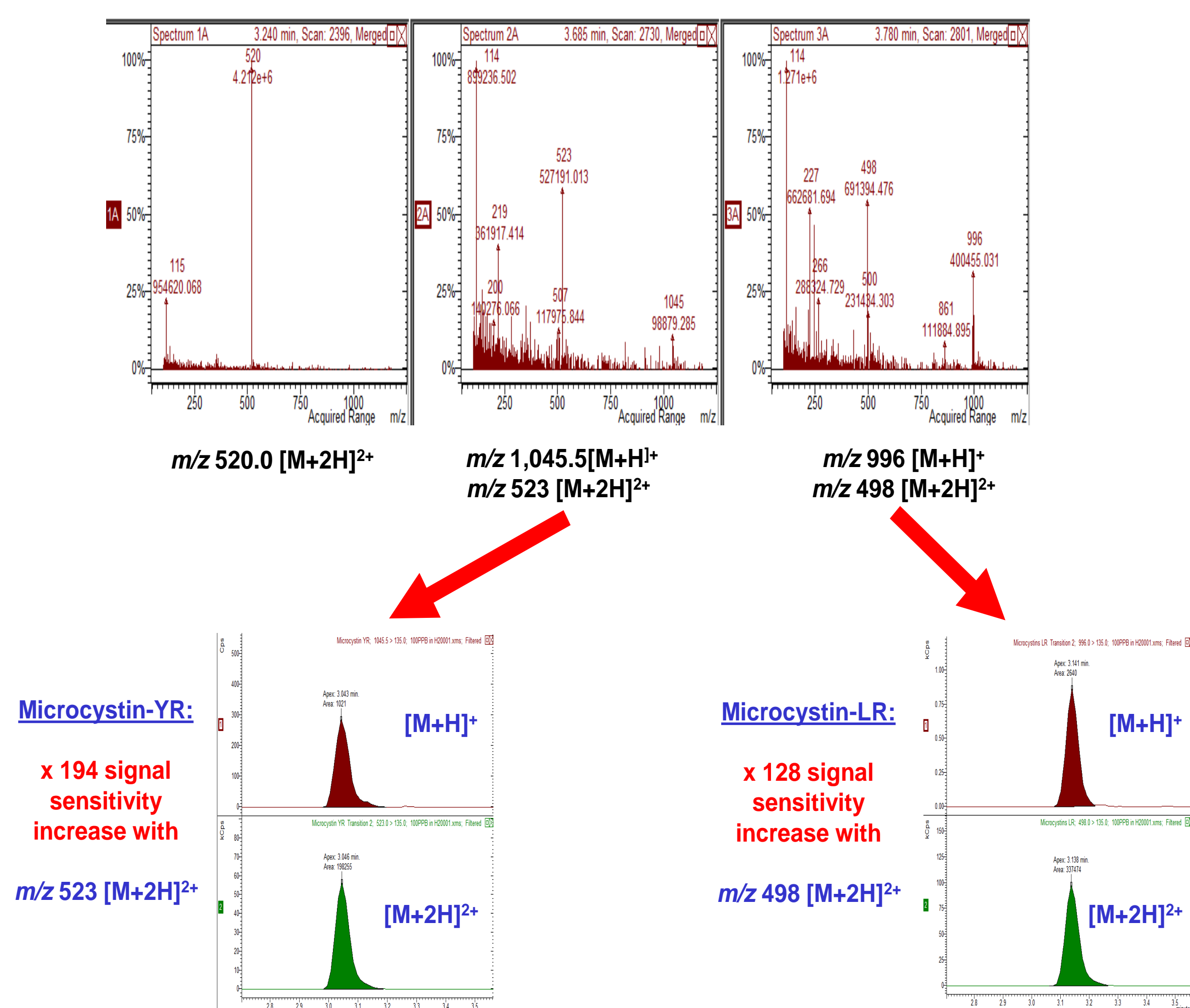

## 5. REPRODUCIBILITY AND CALIBRATION DATA

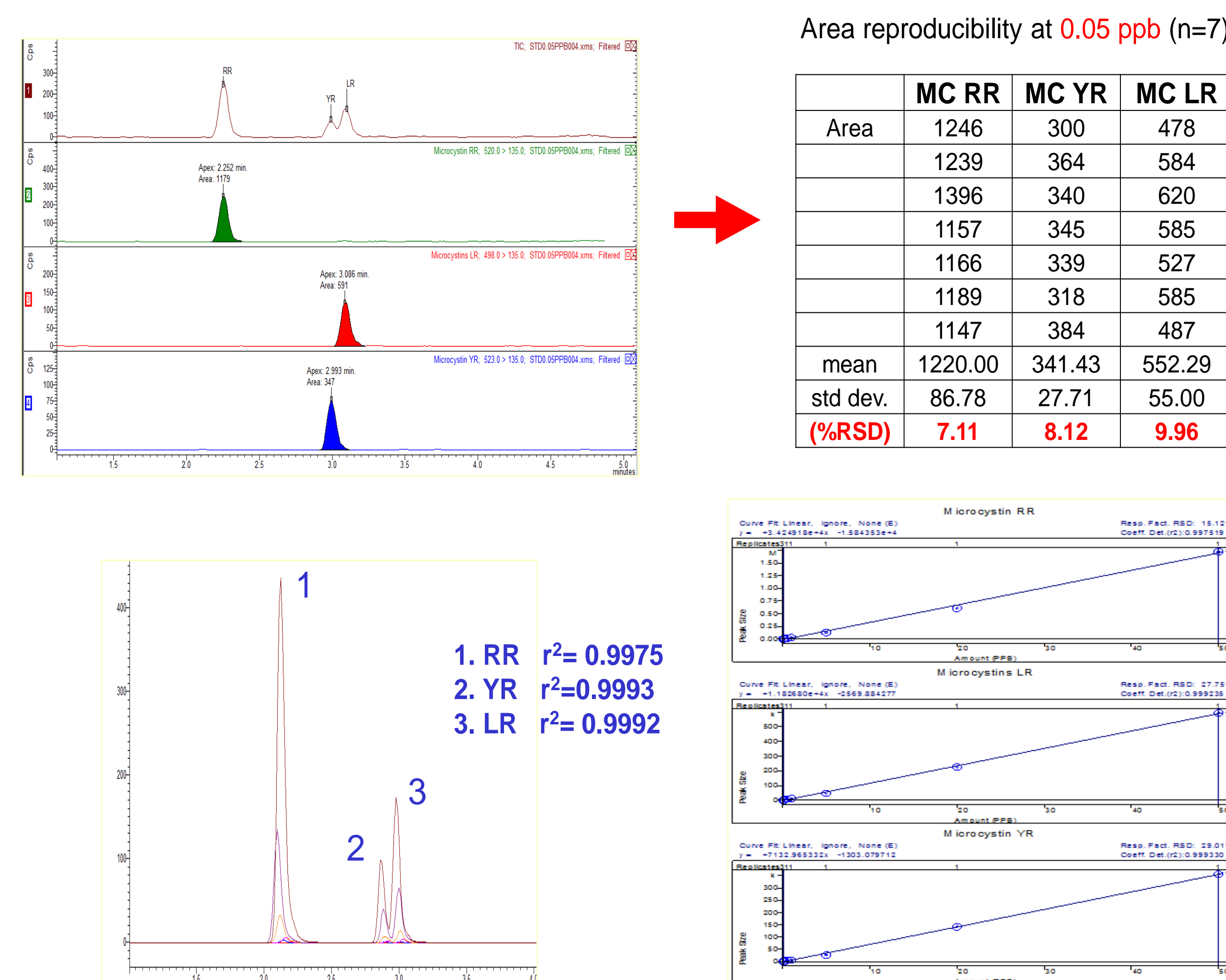

## 3. UHPLC AND MS CONDITIONS

### Bruker Advance™ UHPLC

Column: ACE Excel 2µm C18, 100 x 2.1mm  
Injection volume: 50µL  
Flow rate: 0.4mL/min  
Column temperature: 40°C  
Mobile phase A: 0.1% formic acid (aq)  
Mobile phase B: MeCN

### Gradient conditions:

Time (mins)	%B
0	30
1.0	30
7.0	95
7.1	30
10.0	30

### Bruker EVOQ™

VIP Heated-ESI Temp: 350 °C  
Heated gas: 50 units  
Nebulizer gas: 50 units  
Cone gas temp: 200 °C  
Cone gas: 15 units  
Spray voltage: 4500V (Positive)  
Active exhaust: On  
Collision gas: Argon 1.5 mTorr  
Optimized MRM transitions:

MC RR (MW: 1038) transition: 520 → 135 (CE at 24V)  
MC LR (MW: 995) transition: 498 → 135 (CE at 11V)  
MC YR (MW: 1045) transition: 523 → 135 (CE at 9V)



## 6. SUMMARY AND CONCLUSIONS

- A new method was established using an ACE Excel 2µm C18 column and Bruker Advance / EVOQ UHPLC-MS/MS instrumentation.
- The method was found to be reproducible and sensitive at very low levels (0.05ppb) for microcystins in water.
- Rapid separation and elution of Microcystin-LR, Microcystin-RR and Microcystin-YR was achieved with an excellent linear range of 0.05ppb – 50ppb.
- MS optimization work indicated significantly enhanced sensitivity was possible using doubly charged ions.

ACE® is a registered trademark of Advanced Chromatography Technologies Limited. Excel™ is a trademark of Advanced Chromatography Technologies Limited. Advanced Chromatography Technologies Limited acknowledges the registered and unregistered trademarks of Bruker and has no affiliation with the company.

All registered and unregistered trademarks are recognised. All data reproduced with permission of Bruker from an original slide deck by Kyung Hoon Cha, Bruker Korea. Contact Bruker for more information.