



Mycotoxins / Aflatoxins from Peppers

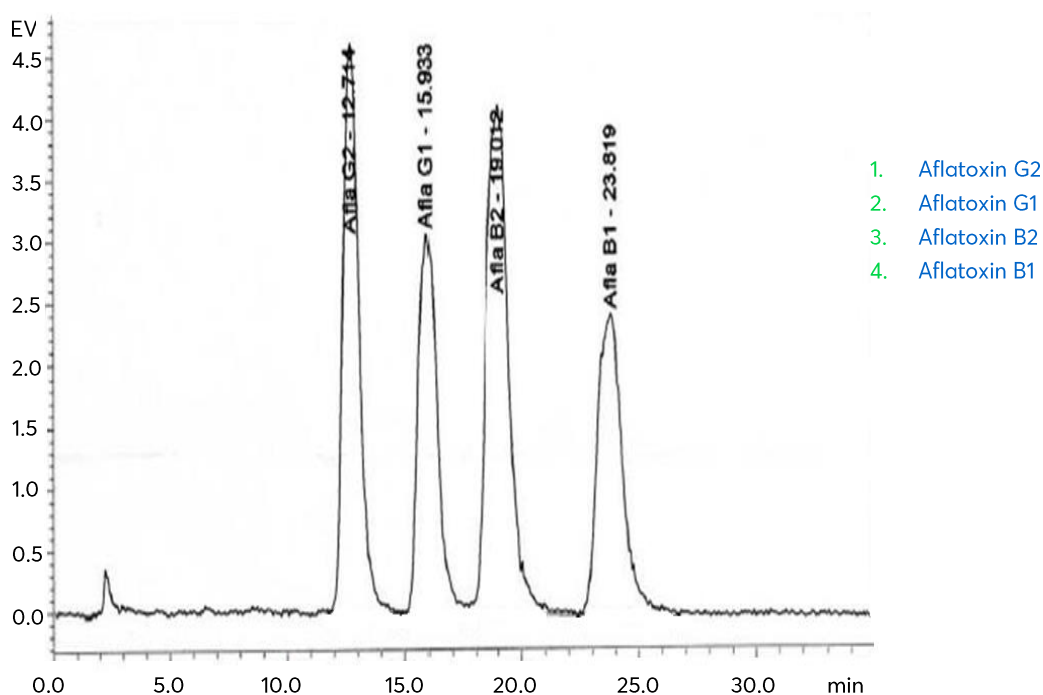


Method Details

CONDITIONS


Column:	Avantor® ACE® 3 C18-PFP
Particle Size:	3 µm
Dimensions:	150 x 4.6 mm
Mobile Phase:	H ₂ O/MeOH (60:40 v/v)
Flow Rate:	1 mL/min
Injection:	100 µL
Temperature:	45 °C
Detection:	Fluorescence, Ex λ 362 nm, Em λ 425 nm

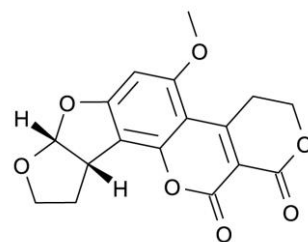


Aflatoxin G2

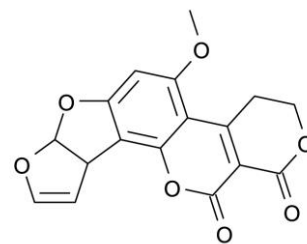

Aflatoxin G1

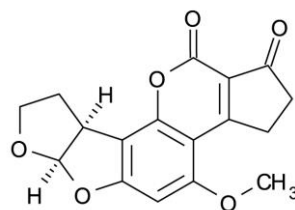

Aflatoxin B2

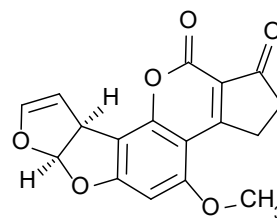

Aflatoxin B1

ORDERING TABLE

Product	Details	Size	Part Number
Avantor® ACE® 3 C18-PFP	HPLC Column	150 x 4.6 mm	ACE-1110-1546

Reproduced with permission of Public Health Laboratory of Soria, Spain