


Nitrosamine Contaminants in Valsartan API

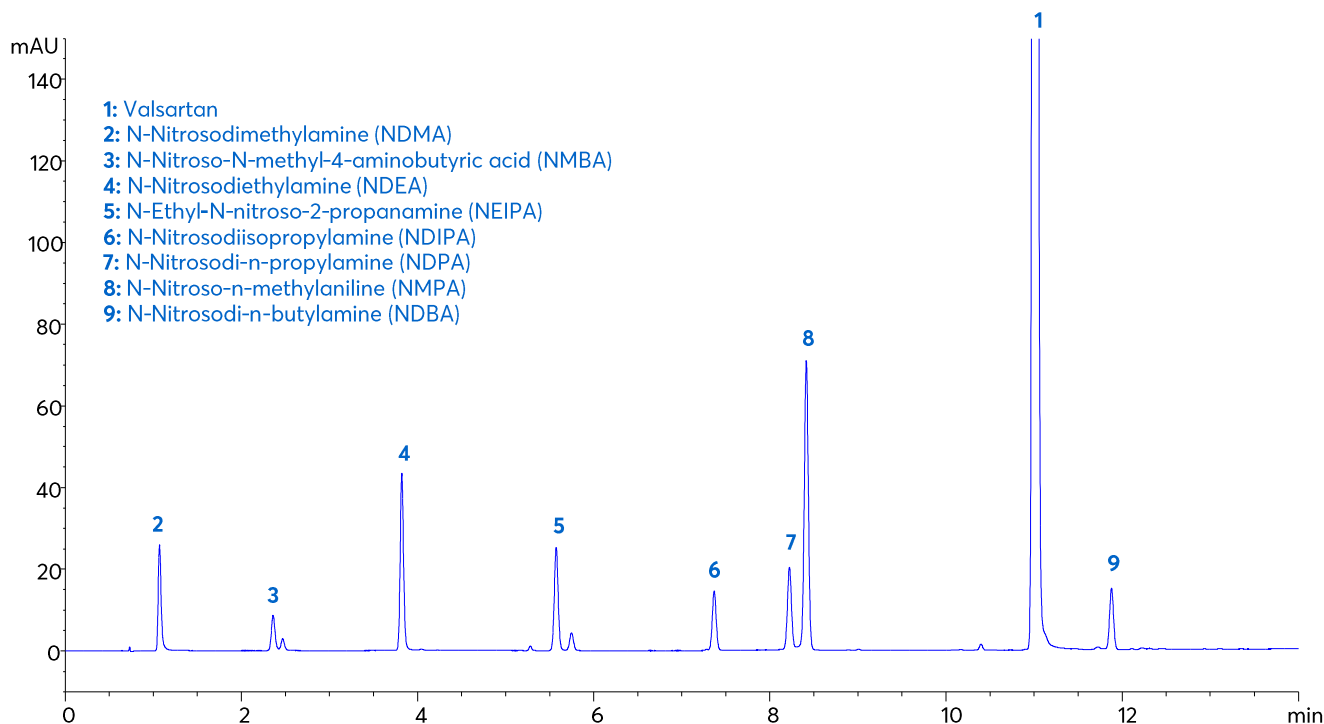

Method Details

CONDITIONS

Column: Avantor® ACE® UltraCore C18

Dimensions: 100 x 3.0 mm

Mobile Phases: A: 20 mM KH₂PO₄ pH 2.7 in H₂O
B: 20 mM KH₂PO₄ pH 2.7 in MeCN/H₂O 70:30 v/v

Gradient:

Time (mins)	%B
0	4
1	4
15	95
17	95
17.5	4
25	4

Flow Rate: 0.6 mL/min

Injection: 1 µL

Temperature: 20 °C

Detection: UV, 254 nm

System Dwell Volume: 204 µL

ORDERING TABLE

Product	Details	Size	Part number
Avantor® ACE® UltraCore C18	HPLC column	100 x 3.0 mm	CORE-35F-1030
Acetonitrile	VWR CHROMANORM® gradient grade	2.5 L	83639.320
Water	VWR CHROMANORM® gradient grade	2.5 L	83650.320
Potassium dihydrogen phosphate	VWR HiPerSolv CHROMANORM® for HPLC	500 g	153184U