


Nucleic Acids / Disease Biomarker Profiling I

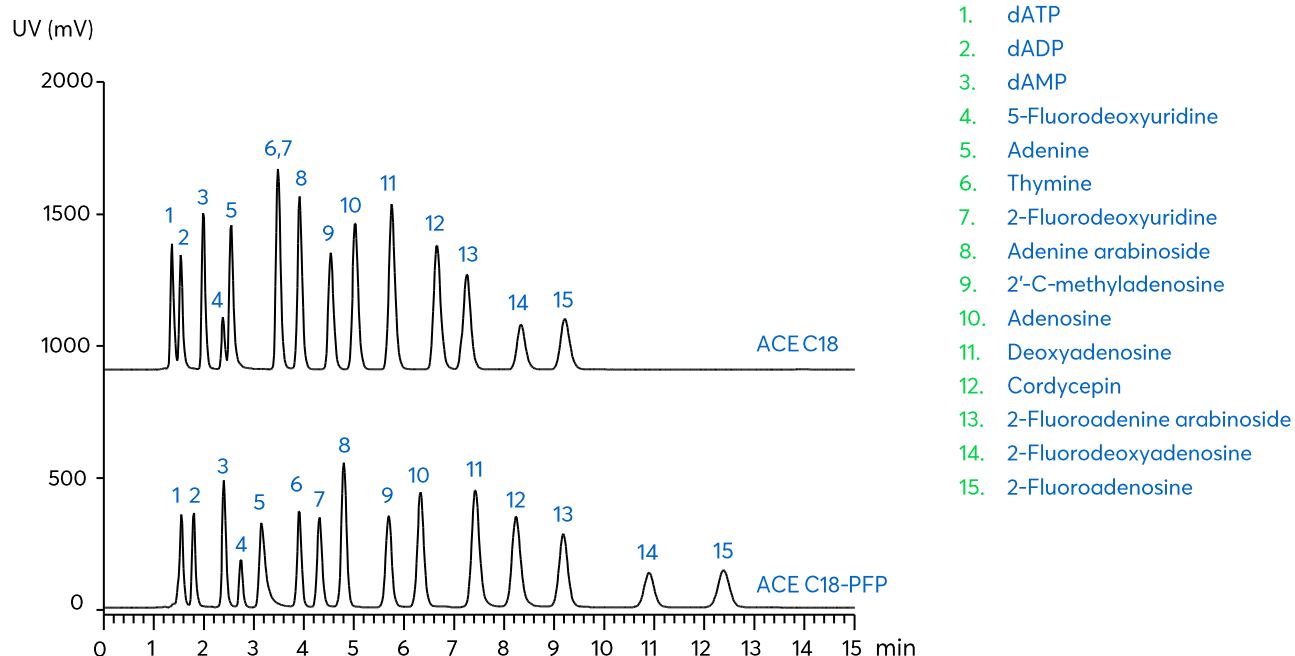

Method Details

CONDITIONS

Column:	Avantor® ACE® 3 C18-PFP Avantor® ACE® 3 C18
Particle Size:	3 µm
Dimensions:	100 x 4.6 mm
Mobile Phase:	33 mM potassium phosphate pH 6.2 with KOH/MeOH (88:12 v/v)
Flow Rate:	1 mL/min
Temperature:	Ambient
Detection:	UV, 260 nm

ORDERING TABLE

Product	Details	Size	Part Number
Avantor® ACE® 3 C18-PFP	HPLC Column	100 x 4.6 mm	ACE-1110-1046
Avantor® ACE® 3 C18	HPLC Column	100 x 4.6 mm	ACE-111-1046

Reproduced with permission of Department of Medical Biochemistry and Biophysics, Umeå University, Sweden