

### Conditions

Column: ACE 5 C18-AR  
 Dimensions: 150 x 4.6 mm  
 Part Number: ACE-129-1546  
 Mobile Phase: A: 5 mM formic acid in H<sub>2</sub>O  
 B: 5 mM formic acid in MeOH

Time (mins)	%B
0.0	3
5.0	100
5.5	100
6.0	3
Post time 2.5 minutes	

Flow Rate: 1.5 mL/min  
 Injection: 5 µL  
 Temperature: 40 °C  
 Detection: UV, 254 nm

1. Paracetamol
2. Hydrochlorothiazide
3. Methylphenylsulfoxide
4. Methylphenylsulfone
5. Aspirin
6. Phenacetin
7. 1,3-Dinitrobenzene
8. 1,2,4-Trimethoxybenzene
9. Ethylbenzoate
10. Nimesulide
11. Ibuprofen
12. Indomethacin
13. Mefenamic acid

