


Application note #2290

Phenol and Phenoxy Acid Herbicides

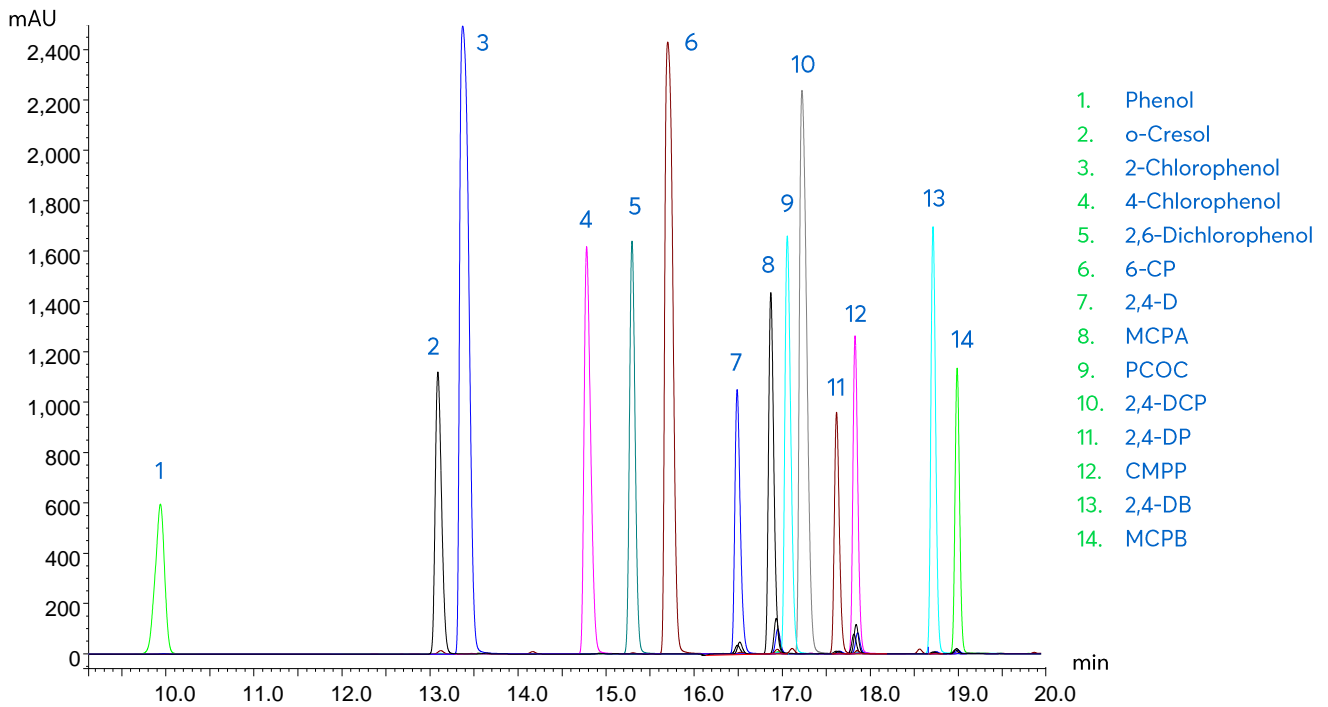

Method Details

CONDITIONS

Column: Avantor® ACE® 3 C18-PFP
 Particle Size: 3 µm
 Dimensions: 150 x 4.6 mm
 Mobile Phase: A: 0.1% formic acid in H₂O
 B: MeOH

Time (mins)	% B
0.0	10
20.0	100

Flow Rate: 1 mL/min
 Injection: 10 µL
 Temperature: 35 °C
 Detection: UV, 280 nm

ORDERING TABLE

Product	Details	Size	Part Number
Avantor® ACE® 3 C18-PFP	HPLC Column	150 x 4.6 mm	ACE-1110-1546

Reproduced with permission of Nufarm UK Ltd