

# Sennosides in Traditional Chinese Medicine

**ACE**<sup>®</sup>  
Ultra-inert  
UHPLC & HPLC Columns



Application #AN1390

## Conditions


Column: ACE 3 C18-PFP  
Dimensions: 150 x 4.6 mm  
Part Number: ACE-1110-1546  
Mobile Phase: A: 0.75% acetic acid in H<sub>2</sub>O  
B: MeCN/MeOH (90:10 v/v)

Time (mins)	%B
0	9
23	28
40	28

Flow Rate: 0.6 mL/min  
Temperature: 35 °C  
Detection: UV, 271 nm  
Sample: Herbal tea bag containing Folium Sennae, Peppermint, Folium Mori, Folium Nelumbinis, Glycyrrhiza Uralensis and Lalang Grass Rhizome

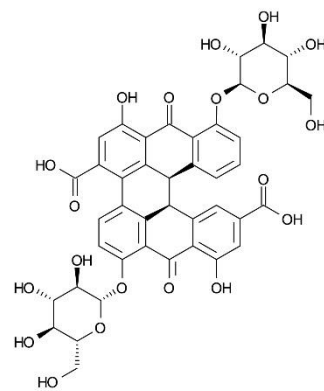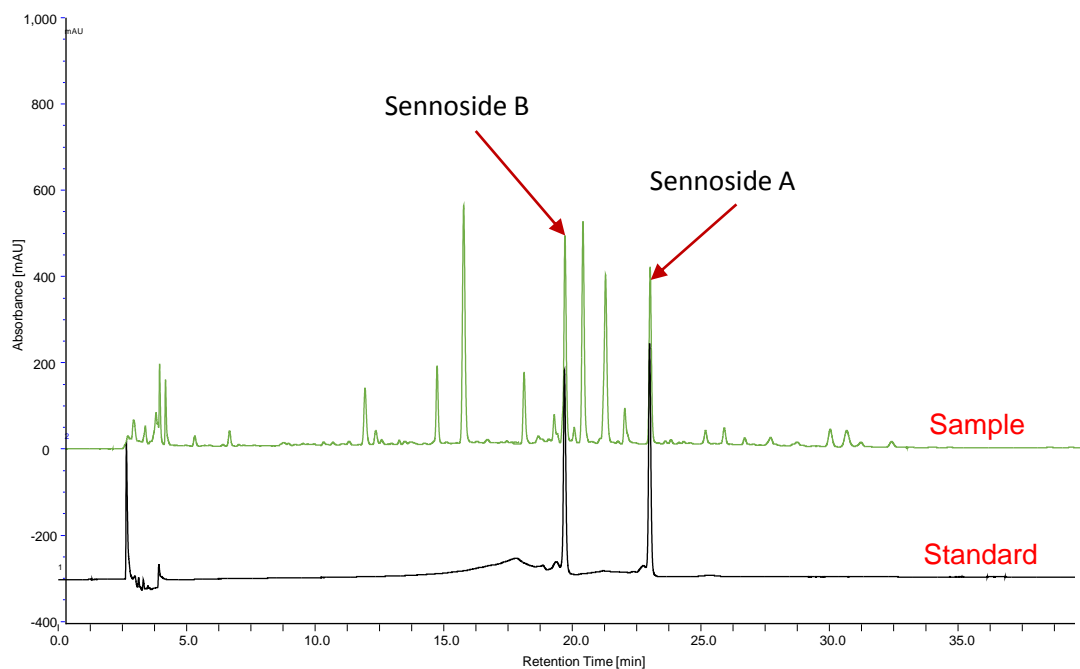

Sennoside A

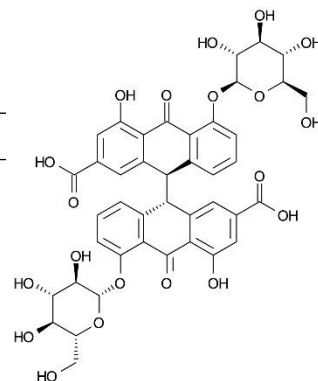

Sennoside B

Reproduced with permission of Vercotech Inc., Taiwan

© 2015 ACT Ltd. All rights reserved. All products are available worldwide.

For more information contact your local ACE distributor or visit  
[www.ace-hplc.com](http://www.ace-hplc.com) or email: [info@ace-hplc.com](mailto:info@ace-hplc.com)

**ACE**<sup>®</sup>  
UHPLC & HPLC  
Columns