

Steroids Separation using Enhanced Polar Selectivity

ACE[®]
Ultra-inert
UHPLC & HPLC Columns

Application #AN2470

Conditions

Column: ACE Excel 3 C18-Amide
Dimensions: 50 x 2.1 mm
Part Number: EXL-1112-0502U
Mobile Phase: A: 0.1% formic acid in H₂O
B: 0.1% formic acid in MeCN
Time (mins) %B
0 25
10 80
Flow Rate: 0.4 mL/min
Temperature: 20 °C
Detection: UV, 260 nm

