



Application #AN3680

Conditions


Column:	ACE 5 C18
Dimensions:	150 x 3.0 mm
Part Number:	ACE-121-1503
Mobile Phase:	10 mM oxalic acid pH 2.9/MeCN (80:20 v/v)
Flow Rate:	0.5 mL/min
Injection:	20 µL
Temperature:	Ambient
Detection:	UV-Vis, 350 nm

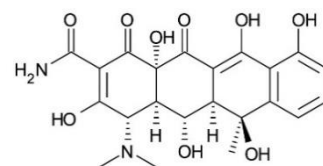


1. Oxytetracycline

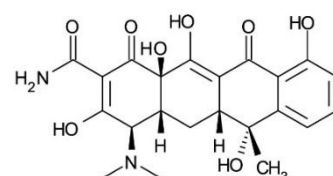


2. Tetracycline

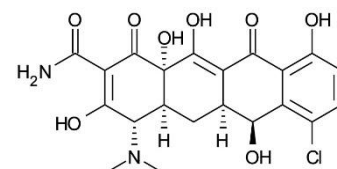

3. Demeclocycline

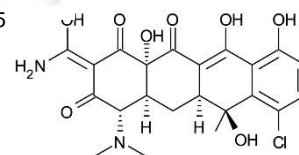

4. Chlortetracycline

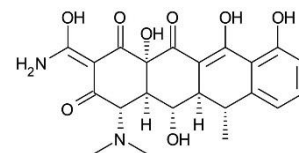

5. Doxycycline

