


Conditions


Column: ACE Excel 2 SuperC18
Dimensions: 100 x 3.0 mm
Part Number: EXL-1011-1003U
Mobile Phase: A 20 mM ammonium formate pH 3.0 in H₂O
B 20 mM ammonium formate pH 3.0 in MeOH/H₂O (9:1 v/v)

Time (mins)	%B
0.0	50
6.0	70
7.0	70
7.5	50


Flow Rate: 1.2 mL/min
Injection: 2 µL
Temperature: 40 °C
Detection: UV, 260 nm

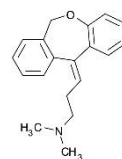


1. Doxepin + isomer

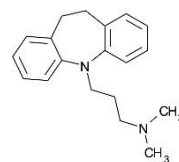


2. Imipramine

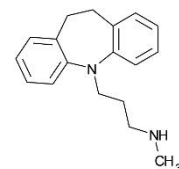


3. Desipramine

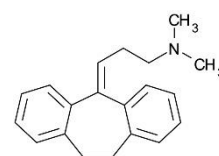

4. Amitriptyline

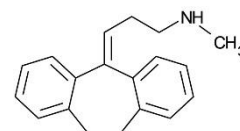

5. Nortriptyline

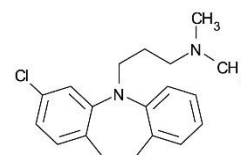

6. Clomipramine

