

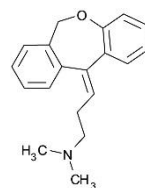


## Conditions


Column:	ACE Excel 2 C18-PFP
Dimensions:	100 x 3.0 mm
Part Number:	EXL-1010-1003U
Mobile Phase:	20 mM ammonium formate pH 3.0 in MeOH/H <sub>2</sub> O (54:46 v/v)
Flow Rate:	1.2 mL/min
Injection:	2 µL
Temperature:	40 °C
Detection:	UV, 260 nm

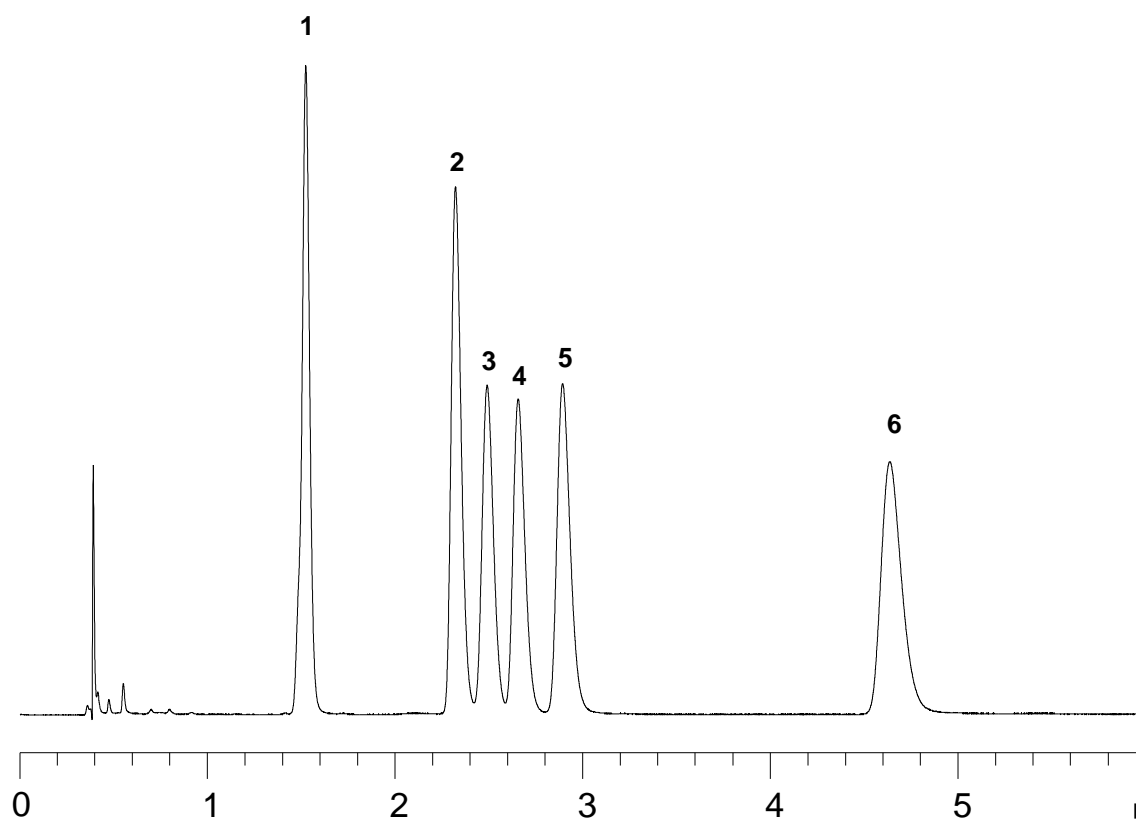


1. Doxepin

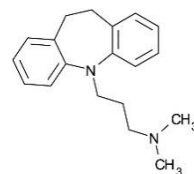


2. Imipramine

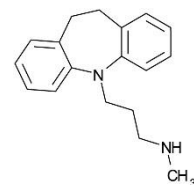


3. Desipramine

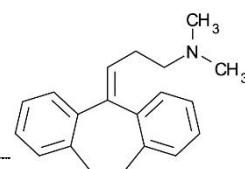

4. Amitriptyline

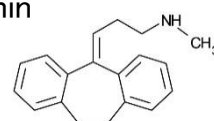

5. Nortriptyline

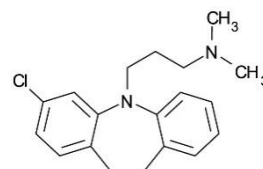

6. Clomipramine