

1. Doxepin
2. Protriptyline
3. Imipramine
4. Nortriptyline
5. Amitriptyline


Conditions (100x3mm):

60/40 MeCN/20mM Phosphate
(pH7.0)


Wavelength: 254nm

Flowrate: 0.4ml/min

Temperature: 25C

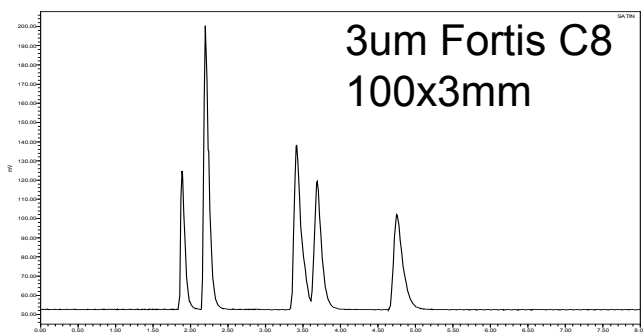

**Selectivity –
Less
Retention**

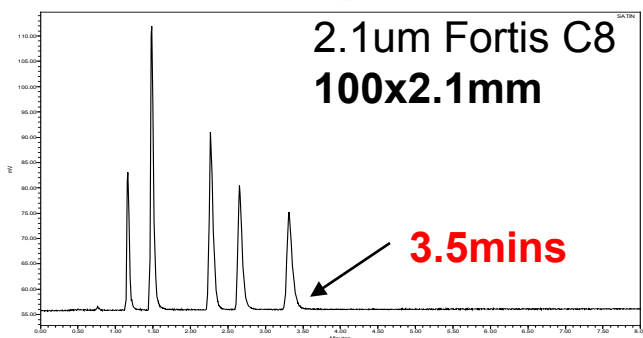

**Efficiency –
More
Resolution**

Conditions (100x2.1mm):

60/40 MeCN/20mM Phosphate
(pH7.0)

Wavelength: 254nm

Flowrate: 0.4ml/min

Temperature: 40C