
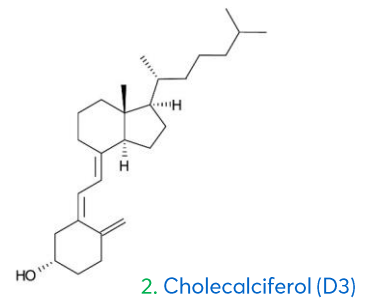
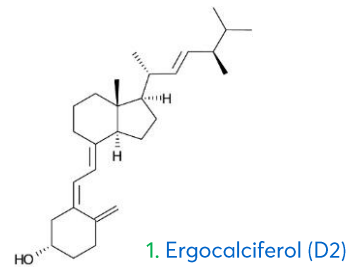


Vitamin D2/D3

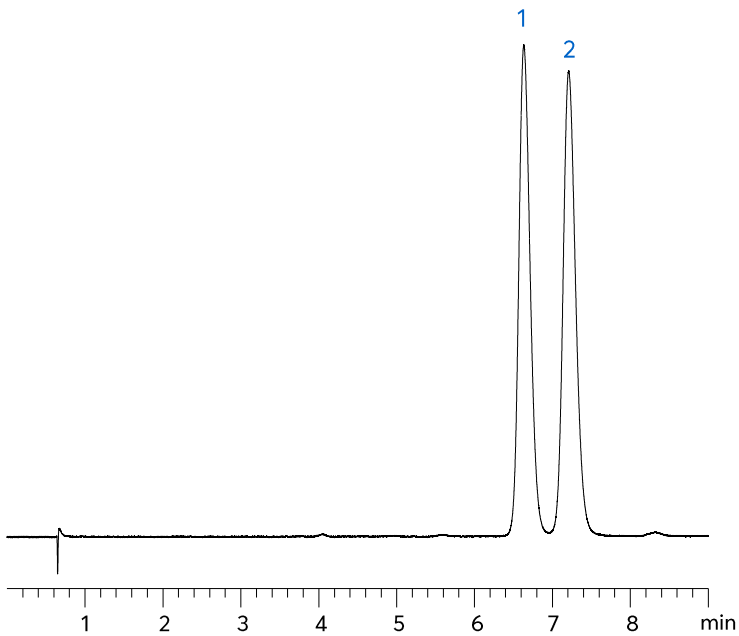

Method Details

CONDITIONS

Column:	Avantor® ACE® Excel® 2 C18-Amide
Particle Size:	2 µm
Dimensions:	50 x 3.0 mm
Mobile Phase:	100% MeCN
Flow Rate:	0.43 mL/min
Injection:	2 µL
Temperature:	20 °C
Detection:	UV, 265 nm

ORDERING TABLE

Product	Details	Size	Part Number
Avantor® ACE® Excel® 2 C18-Amide	UHPLC Column	50 x 3.0 mm	EXL-1012-0503U
Acetonitrile	VWR CHROMANORM® super gradient grade	2.5 L	83639.320